# IKA® Vortex Genius 3

Betriebsanleitung Ursprungssprache	DE	03
Operating instructions	EN	07
Mode d'emploi	FR	11
Руководство пользователя	RU	15

Indicaciones de seguridad	ES	<b>19</b>	Biztonsági utasítások	HU	<b>29</b>
Veiligheidsinstructies	NL	<b>20</b>	Varnostna opozorila	SL	<b>30</b>
Norme di sicurezza	IT	<b>21</b>	Bezpečnostné pokyny	SK	<b>31</b>
Säkerhetsanvisningar	SV	<b>22</b>	Ohutusjuhised	ET	<b>32</b>
Sikkerhedshenvisninger	DA	<b>23</b>	Drošības norādes	LV	<b>33</b>
Sikkerhedsanvisninger	NO	<b>24</b>	Nurodymai dėl saugumo	LT	<b>34</b>
Turvallisuusohjeet	FI	<b>25</b>	Инструкции за безопасност	BG	<b>35</b>
Instruções de serviço	PT	<b>26</b>	Indicații de siguranță	RO	<b>36</b>
Wskazówki bezpieczeństwa	PL	<b>27</b>	Υποδείξεις ασφάλειας	EL	<b>37</b>
Bezpečnostní pokyny	CS	<b>28</b>			

## Bedienelemente / Control elements / Éléments de réglage / Коммутационные элементы


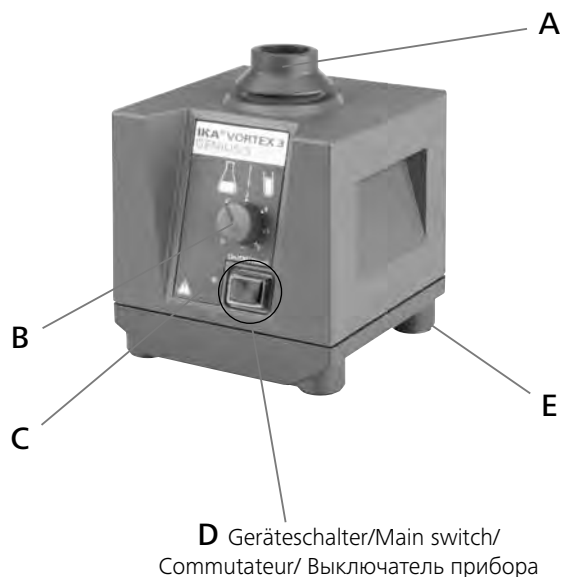
Fig. 1

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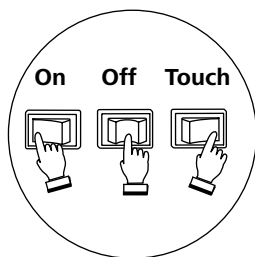	Page		
Control elements	02	Correct use	08
EC - Declaration of Conformity	02	Operation	08
Contents	07	Accessories	09
Explication of warning symbols	07	Maintenance	10
Safety instructions	07	Warranty	10
Unpacking	08	Technical data	10

## Explication of warning symbols

General hazard.

This symbol identifies information **that is of vital importance for protecting your health and safety**. Disregarding this information may lead to health impairment and injuries.

This symbol identifies information **that is of importance for the technically correct functioning of the system**.

Disregarding this information can result in damage to the instrument or to system components.

## Safety instructions

### General information

- **Read the operation instructions completely before starting up and follow the safety instructions.**
- Keep the operation instructions in a place where they can be accessed by everyone.
- Ensure that only trained staff work with the appliance.
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- Socket must be earthed (protective ground contact).

### Device design

- 
  - **Do not** operate the appliance in explosive atmospheres, with hazardous substances or under water.
- Set up the appliance in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.
- The feet of the appliance must be clean and undamaged.
- The voltage stated on the type plate must correspond to the mains voltage.
- The socket for the mains cord must be easily accessible
- Check the appliance and accessories for damage before each use. Do not use damaged components..
- Firmly secure the accessories and vessels in place, otherwise shaking vessels could be damaged or projected out
- Position one single shaking vessel in the centre and several shaking vessels so that they are evenly spread out.
- Sharp edged vessels cause wear debris on the attachments.

### Permissible medium / contaminants / side reactions

- 
  - Beware of hazards due to:
    - flammable materials
    - glass breakage

- Only process media that will not react dangerously to the extra energy produced through processing. This also applies to any extra energy produced in other ways, e.g. through light irradiation.
- The appliance may heat up when in use.

### Power supply / Switching off

- The appliance starts up again automatically following a cut in the power supply.

### Experimental procedures

- Wear your personal protective equipment in accordance with the hazard category of the medium to be processed. Otherwise there is a risk of:
  - splashing liquids
  - projectile parts
  - body parts, hair, clothing and jewellery getting caught.
- Position the knob (B) at the left stop before starting up. Gradually increase the speed.

- Keep hand/arm vibration to a minimum (Directive 2002/44/EC) when operated manually:
  - Keep the sample container as vertical as possible.
  - Only exert the minimum pressure necessary for the mixing process.
- If the shaking movement of the appliance increases (resonance), reduce the speed or pass through the critical phase as quickly as possible.

- Reduce the speed if
  - the medium splashes out of the vessel because the speed is too high
  - the appliance is not running smoothly
  - the appliance begins to move around because of dynamic forces.

#### Accessories

- Safe operation is only guaranteed with the accessories described in the "Accessories" chapter.

- Always disconnect the plug before attaching accessories.
- Accessories must be securely attached to the device and cannot come off by themselves. The centre of gravity of the assembly must lie within the surface on which it is set up.

#### For protection of the equipment

- The appliance may only be opened by experts.
- Protect the appliance and accessories from bumps and impacts.

## Unpacking

### • Unpacking

- Please unpack the device carefully
- In the case of any damage a report must be sent immediately (post, rail or forwarder).

### • Scope of delivery

- Shaker **Vortex Genius 3**
- Standard attachment VG 3.1
- Operating instruction

## Correct use

### • Use

- For mixing liquids
  - Touch mode for single test tubes
  - Continuous mode for single or several vessels

### • Range of use (only indoors)

- Laboratories      - Schools
- Pharmacies

This device is suitable for use in all areas except:

- Residential areas
- Areas that are connected directly to a low-voltage supply network that also supplies residential areas.

The safety of the user cannot be guaranteed if: the appliance is operated with accessories that are not supplied or recommended by the manufacturer, the appliance is operated improperly according to the manufacturer's specifications or the appliance or printed circuit board are modified by third parties.

## Commissioning

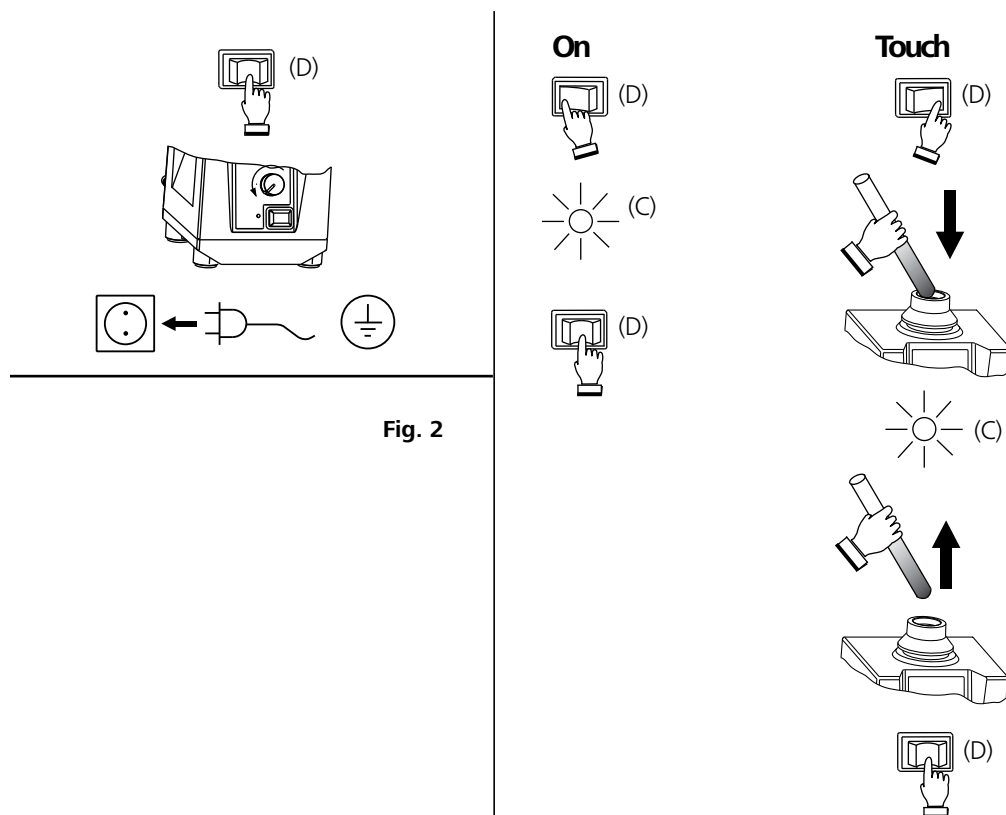
Fig. 2

## Accessories

Possible uses and permitted speed ranges of attachments:

Designation	Description	Id. - No.	Touch Mode	Continuous Mode	Range of speed
 VG 3.1 Standard attachment	For test tubes and small vessels upto diameter 30 mm	33 412 00	X	X	
 VG 3.2 One-hand attachment	88 mm-plate attachment with rubber pad	33 423 00	X	X	
 VG 3.3 150 mm-Universal attachment	150 mm-plate attachment with rubber pad	33 424 00	-	X	
 VG 3.31 Test tube insert	For 54 Eppendorf tubes for VG 3.3	33 443 00	-	X	
 VG 3.32 Test tube insert	For 18 test tubes 10 mm, for VG 3.3	33 439 00	-	X	
 VG 3.33 Test tube insert	For 12 test tubes 12 mm, für VG 3.3	33 440 00	-	X	
 VG 3.34 Test tube insert	For 8 test tubes 16 mm, for VG 3.3	33 441 00	-	X	
 VG 3.35 Test tube insert	For 8 test tubes 20 mm, for VG 3.3	33 442 00	-	X	
 VG 3.36 Erlenmeyer flask attachment	For 1 Erlenmeyer-/ round-bottomed flask 100-250 ml, for VG 3.3	33 421 00	-	X	
 VG 3.37 Microtiter plate insert-	For 1 microtiter plate, for VG 3.3	33 444 00	-	X	

### Changing attachments

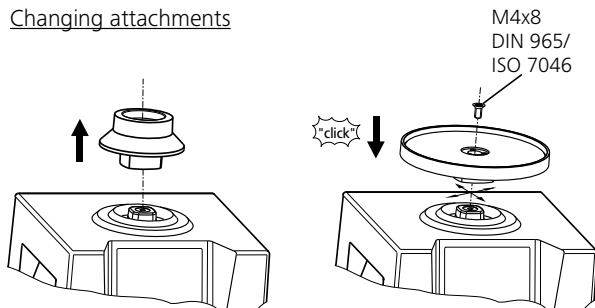
Fig. 3

### Using inserts


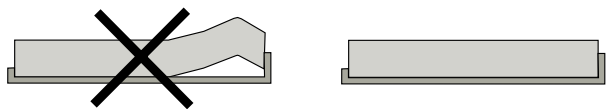
Fig. 4

## Maintenance

The device is maintenance-free. It is subject only to the natural wear and tear of components and their statistical failure rate.

### Cleaning

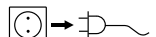For cleaning, disconnect the main plug!

Only use cleaning agents recommended by **IKA®**.

### Contamination Cleaning agents

Dyes	isopropyl alcohol
Building materials	water containing tenside/isopropyl alcohol
Cosmetics	water containing tenside/isopropyl alcohol
Foodstuffs	water containing tenside
Fuels	water containing tenside

- When cleaning, please do not allow moisture to get into the unit.
- Wear the proper protective gloves during cleaning of the devices.
- Before performing a non-recommended method of cleaning or decontamination, the user must ascertain with **IKA®** that this method does not destroy the instrument.

### Spare parts order

When ordering spare parts, please give:

- Machine type
- Serial number, see type plate
- Software Version (second displayed value when the device is on)
- Item and designation of the spare part please see

### Repair

**Please send instrument in for repair only after it has been cleaned and is free from any materials which may constitute a health hazard.**

For this you should request the “**Decontamination Clearance Certificate**” from **IKA®**, or use the download printout of it from the **IKA®** website

Return the instrument in its original packaging. Storage packaging is not sufficient. Also, please use suitable shipping package materials.

## Warranty

In accordance with **IKA®** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine directly to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

## Technical data

Operating voltage	<b>VAC</b>	220 – 240 ± 10%
Design frequency	<b>Hz</b>	50/60
Nominal voltage	<b>V/Hz</b>	230/50
Input power	<b>W</b>	60
Output power	<b>W</b>	10
Speed	<b>rpm</b>	0/ 500 – 2500 infinitely adjustable
Speed adjustment		analog (infinitely) rotating knob in front of device
Speed display		Scale 0 – 6
Drive		Shadepole motor
Agitation stroke	<b>mm</b>	4
Shaking motion		horizontal, approximately circular
Perm. duration of operation	<b>%</b>	100
Perm. ambient temperature	<b>°C</b>	+5 to +40
Perm. relative humidity	<b>%</b>	80
Protection type acc. to DIN EN 60529		IP 21
Protection class		I (protection earth)
Overvoltage category		II
Contamination level		2
Operation at terrestrial altitude	<b>m</b>	max. 2000
Dimensions	<b>mm</b>	127 x 149 x 136
total W x D x H		
Weight without supported load	<b>kg</b>	4,5
Max. supported load incl. attachment	<b>kg</b>	0,5

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