

# INSTRUCTIONS

---

# MVX10

## Research high-class stereo microscope

This instruction manual is for the EVIDENT MVX Research Macro Zoom System Microscope. To ensure the safety, obtain optimum performance and familiarize yourself fully with the use of this attachment, we recommend you study this manual thoroughly before operating the microscope. Retain this instruction manual in an easily accessible place near the work desk for future reference. For details of products included in the configuration of this system, see "1 Nomenclature" (P.4) and "9 Operation of other modules" (P.24).

Optical Microscope and Accessory

# Contents

Correct assembly and adjustments are indispensable for the microscope to manifest its full performance. If you want to assemble the microscope by yourself, refer to "10 Assembly" on page 33.

<b>Important</b> .....	<b>1</b>
<b>1 Nomenclature</b> .....	<b>4</b>
<b>2 Controls</b> .....	<b>5</b>
3-1 Preparation.....	8
3-2 Observation procedure .....	8
<b>4 Operation</b> .....	<b>10</b>
4-1 Base .....	10
4-2 Microscope body and focusing assembly.....	10
<b>1</b> Adjusting the focus.....	10
<b>2</b> Adjusting the rotation tension of the coarse focus adjustment knob .....	10
<b>3</b> Engaging and disengaging the zooming knob click stop position.....	11
<b>4</b> Adjusting the aperture Iris diaphragm.....	11
<b>5</b> Using the objective correction collar.....	12
4-3 Coaxial fluorescence illuminator.....	13
<b>1</b> Selecting the fluorescence mirror unit.....	13
<b>2</b> Turning the burner ON.....	14
<b>3</b> Opening/closing the shutter .....	15
<b>4</b> Using the field Iris diaphragm .....	15
<b>5</b> Switching the filter slider knob .....	15
4-4 Observation tube .....	16
<b>1</b> Adjusting the tilt.....	16
<b>2</b> Adjusting the interpupillar distance.....	16
<b>3</b> Adjusting the diopter (zoom parfocality adjustment).....	17
<b>4</b> Using the eye shades .....	18
<b>5</b> Mounting the eyepiece micrometer disk.....	18
<b>6</b> Selecting the light path.....	18



7	Switching the monaural/stereo view.....	18
<b>5 Camera observation and photomicrography.....</b>		<b>19</b>
1	Selecting the camera adapter magnification .....	19
2	Attaching the camera adapter .....	19
3	Selecting the camera light path.....	19
4	Adjusting the parfocality between observation image and monitor image .....	20
<b>6 Troubleshooting .....</b>		<b>21</b>
<b>7 Specifications.....</b>		<b>22</b>
<b>8 Optical characteristics .....</b>		<b>23</b>
<b>9 Operation of other modules.....</b>		<b>24</b>
9-1	Revolving nosepiece MVX-2RE .....	24
9-2	Magnification changer MVX-CA2X .....	26
9-3	BX stage adapter type 1 SZX-STAD1.....	27
9-4	Stage adapter type 1 SZH-STAD1 .....	28
9-5	BX stage adapter type 2 SZX-STAD2.....	29
9-6	Gliding stage SZH-SG.....	30
9-7	Cup stage SZH-SC.....	31
10-1	Assembly diagram.....	33
<b>10 Assembly.....</b>		<b>33</b>
10-2	Detailed assembly procedures .....	34
10-3	Centration of the mercury (xenon) burner and field Iris diaphragm.....	42
<b>11 Request to issue the Safety Confirmation .....</b>		<b>46</b>






# Important

## Safety precautions

1. After observation of a specimen that involves the risk of infection, be sure to clean the positions that contacted the specimen to prevent infection.
  - To avoid the risk that the specimen drops and splatters, be sure to remove the specimen before moving the microscope.
  - If a specimen is destroyed due to an erroneous operation, immediately take the infection prevention measures.
  - The microscope becomes unstable when its height is increased by a mounted attachment. To prevent the specimen from dropping if the microscope topples down, be sure to take the toppling prevention measures when mounting an attachment.
2. The applicable ultrahigh-pressure mercury burners are the USH-103OL (Ushio Inc.) and HBO103W/2 (OSRAM), both of which are DC mercury burners available from EVIDENT.
3. Ensure that the burners are mounted and that the cords are connected securely.
4. The inside of the lamp housing is very hot while the burner is on and immediately after it is turned off. Do not open the lamp housing in these periods (see page 40).
5. The power supply unit contains high-voltage parts inside. Do not attempt to disassemble it.
6. Always use the power cord provided by EVIDENT. If no power cord is provided, please select the power cord by referring to the section "PROPER SELECTION OF THE POWER SUPPLY CORD" at the end of this instruction manual. If the proper power cord is not used, EVIDENT can no longer warrant the electrical safety performance of the equipment. Before plugging the power cord into the wall outlet, ensure that the main switch of the power supply unit is set to "○" (OFF).
7. Always ensure that the grounding terminal of the power supply unit is properly grounded. If the equipment is not grounded, EVIDENT can no longer warrant the electrical safety performance of the equipment.
8. Before opening the lamp housing for replacing the burner, be sure to set the main switch to "○" (OFF), unplug the lamp housing output connector of the power supply unit and wait at least 10 minutes or until the burner and lamp housing have cooled down.
9. The top of the lamp housing becomes very hot. To avoid the risk of a fire, never block the ventilation of this part. Make sure you leave a space of 10 cm or more around the lamp housing and power supply unit to dissipate heat.
10. The power cord is also used to shut off the power supply unit in case of emergency by unplugging. To facilitate this, locate the power supply unit or wall outlet so that the power cord connector (on the rear of the power supply unit) or the wall outlet is easily accessible in case of emergency.



### Safety symbols

The following symbols are found on the microscope. Study the meaning of the symbols and always use the equipment in the safest possible manner.

Symbol	Symbol explanation
	Indicates the presence of a high voltage (1 kV or more), which should not be touched to prevent an electric shock.
	Indicates that the surface becomes hot, and should not be touched with bare hands.
	Before use, carefully read the instruction manual. Improper handling could result in injury to the user and/or damage to the equipment.
	Indicates that the main switch is ON.
	Indicates that the main switch is OFF


### Caution indications

Caution indications are affixed at parts where special precaution is required when handling and using the microscope. Always heed the cautions.


Caution indication positions	• Lamp housing (U-LH100HG, U-LH100HGAP0)	[High temperature caution]	
	• Power supply unit (U-RFL-T)	[High voltage caution]	

# 1 Getting ready

**TIP** This instruction manual pertains only to the operating procedures of the Research Macro Zoom System. Also read the instruction manuals for the modules used in combination with the microscope system so that you can understand the comprehensive operating procedures of the system.

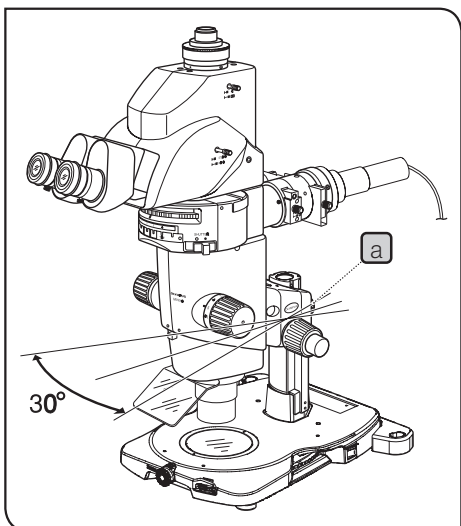

1. A microscope is a precision instrument. Handle it with care and avoid subjecting it to sudden or severe impact. Also be careful in handling the limiting or stopper mechanisms because an excessive force may destroy them.
2. Do not use the microscope where it is subjected to direct sunlight, high temperature and humidity, or vibration. (For the operating environment, see "7 Specifications" (P.22).)
3. To prevent the microscope body from turning over, its pivot angle must be limited to 30° and the tilt angle of the desk surface must be limited to 5° as shown in figure.
4. Care is required when using an auxiliary pillar (SZH-P400/P600) because it increases the height of the microscope body and makes it unstable.
5. To adjust the microscope body height, be sure to hold the focusing assembly with one hand and then loosen the focusing assembly clamping knob **a**. (Use the drop prevention collar (SZX-R) to prevent a hazard from occurring.) Be careful not to pinch your finger during adjustment.
6. Observe the following cautions when operating the coarse focus adjustment knobs or the zooming knobs.

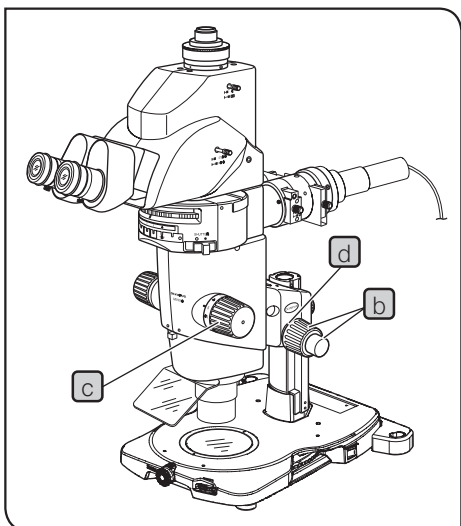

Operation	Manipulated controls	Caution
Focusing	Coarse focus adjustment knobs <b>b</b>	1. the knob hits the upper or lower limiting mechanism violently or it is rotated after it hits a limiting mechanism, the internal mechanism may be damaged.
		2. If the knobs on the left and right are rotated in opposite directions, the internal mechanism will be damaged. (The rotation tension of the knob should be adjusted using the rotation tension adjustment ring <b>d</b> on the knob (see page 10).)
Zooming	Zooming knobs <b>c</b>	1. If the knob hits the upper or lower limiting mechanism violently or it is rotated after it has hit a limiting mechanism, the internal mechanism may be damaged.
		2. If the knobs on the left and right are rotated in opposite directions, the internal mechanism will be damaged.

7. Before moving the microscope, detach the modules including the tilting trinocular head and lamp housing to reduce the total weight. Then hold it by the base, not by the zoom microscope body.

## 2 Maintenance and storage


1. To clean the lenses and other glass components, simply blow dirty away using a commercially available blower and wipe gently using a piece of cleaning paper (or clean gauze).  
If a lens is stained with fingerprints or oil smudges, wipe it gauze slightly moistened with commercially available absolute alcohol




**Since the absolute alcohol is highly flammable, it must be handled carefully.**

**Be sure to keep it away from open flames or potential sources of electrical sparks – for example, electrical equipment that is being switched on or off.**

**Also remember to always use it only in a well-ventilated room.**

2. The equipment uses plastic resins extensively in its external finish. Do not attempt to use organic solvents to clean the non-optical components of the microscope. To clean these components, use a lint-free, soft cloth lightly moistened with a diluted neutral detergent.
3. Never disassemble any part of the microscope as this could result in malfunctions or reduced performance.
4. When not using the microscope, keep it covered with the dust cover provided. Ensure that the lamp housing is cool before covering the microscope.
5. When the hour counter on the power supply unit indicates 300 hours (USH-103OL, HBO103W/2), set the main switch to "O" (OFF) for safety, wait for more than 10 minutes and then replace the burner (see page 40). Unlike the fluorescent lamps, the mercury burner seals high-pressure gas inside. If it used after the specified service life has been exceeded, the glass tube may be distorted accumulatively and may eventually burst, though this happens very rarely. The used mercury burner should be disposed of as an industrial waste. If you cannot dispose of it properly, contact EVIDENT.
6. When disposing of the microscope. Check the regulations and rules of your local government and be sure to observe them. To dispose of only the gas spring  used in the counterbalance of the focusing assembly, follow the precautions provided with the gas spring.

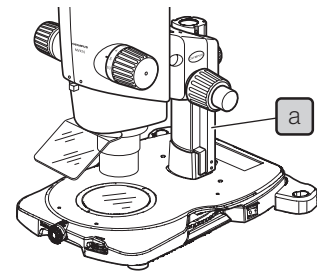

## 3 Caution

If the microscope is used in a manner not specified by this manual, the safety of the user may be imperiled. In addition, the microscope may also be damaged. Always use the microscope as outlined in this instruction manual.

The following symbols are used to set off text in this instruction manual.



**CAUTION** : Indicates that failure to follow the instructions in the warning could result in bodily harm to the user and/or damage to equipment (including objects in the vicinity of the equipment).



**NOTE** : Indicates that failure to follow the instructions could result in damage to equipment.



**TIP** : Indicates commentary (for ease of operation and maintenance.)

## 4 Intended use


This instrument has been designed to be used to observe magnified images of specimens in routine and research applications. Do not use this instrument for any purpose other than its intended use.

# 1 Nomenclature

TIP

The module names listed below are typical ones that can be used with the system, and the illustration shows the system composed of modules marked ●. Certain other modules are also usable even when they are not mentioned below. For these modules, refer to the latest catalogues or contact EVIDENT.


For the modules marked\*, refer to their instruction manuals.

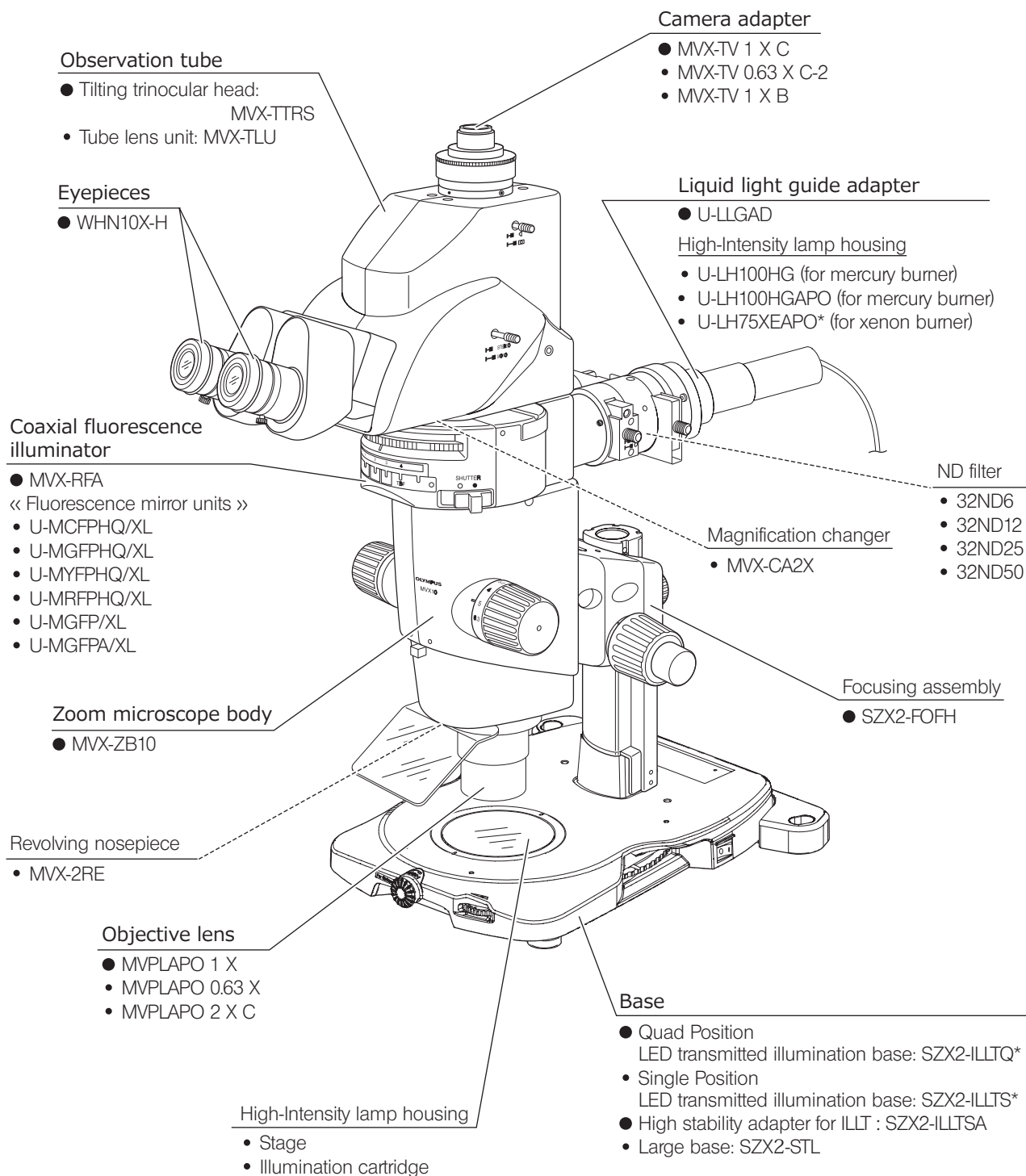

EVIDENT classifies MVX-ZB10 as an optical microscope and other units as optical microscope accessories.

# 2 Controls


**TIP** If the microscope is not yet assembled, go to “10 Assembly” on page 33. The following illustration shows the systems with the modules enclosed in    For other ancillary modules, see subsequent pages. “10 Assembly” (P.33).

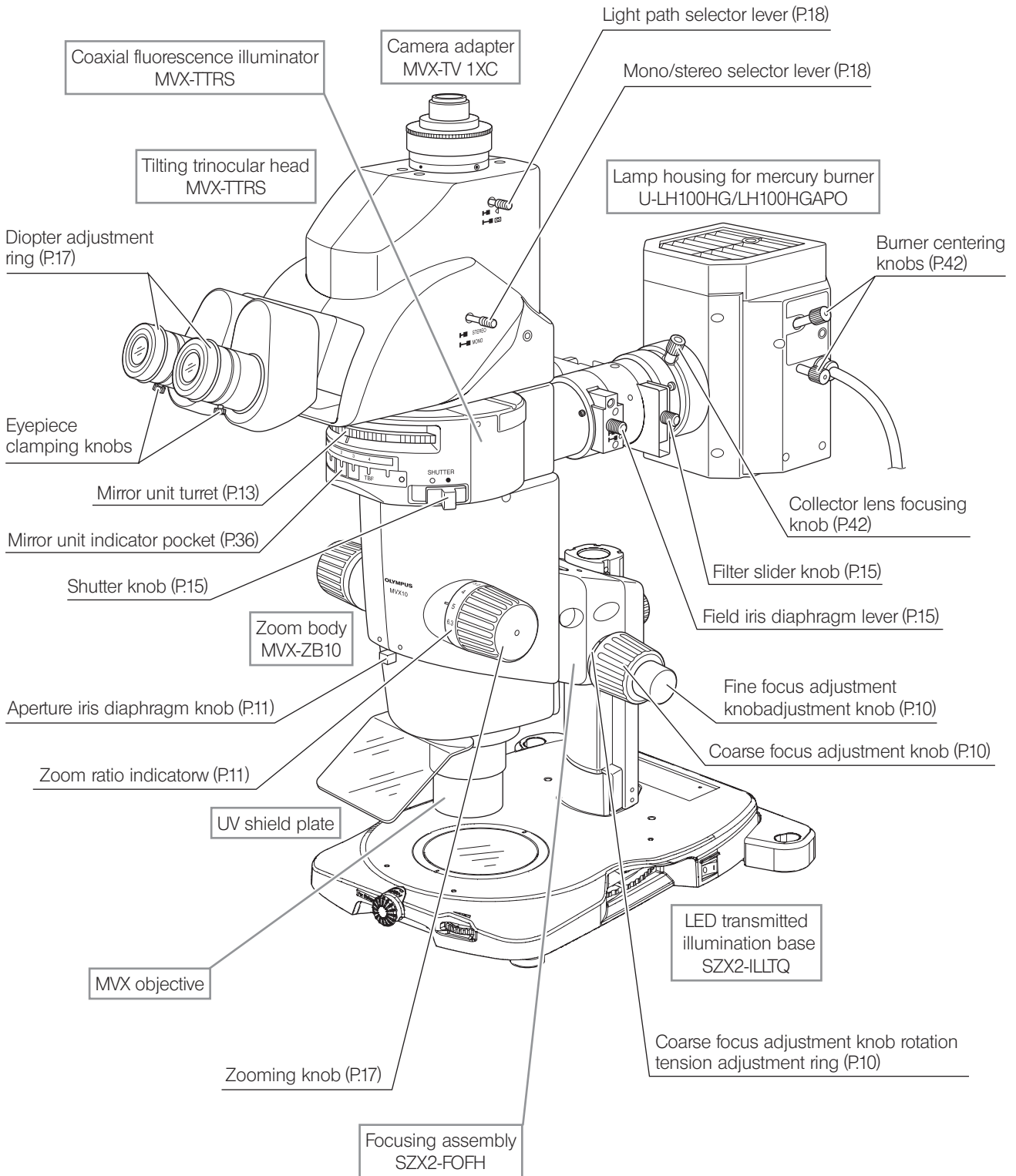

Fluorescence mirror units  
 U-MCFPHQ/XL  
 U-MGFPHQ/XL  
 U-MYFPHQ/XL  
 U-MRFPHQ/XL  
 U-MGFP/XL  
 U-MGFPA/XL


- TIP**
- Up to three fluorescence mirror units can be mounted by placing them in every other dovetail of the six dovetails of the MVX-RFA.
  - When fabricating an original fluorescence unit by your own, it is recommended to use the U-MF/XL mirror unit frame (which does not incorporate a filter). (Page 37) Use the blank indicator sheets provided with the illuminator to write the original mirror unit name.

**NOTE** Each fluorescence mirror unit is composed of the dichroic mirror, barrier filter and excitation filter that match a specific fluorescent protein.

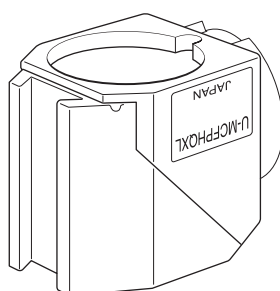

LED and LDP light source  
 U-LGPS

**TIP** For details, see the instruction manual provided with the U-LGPS.

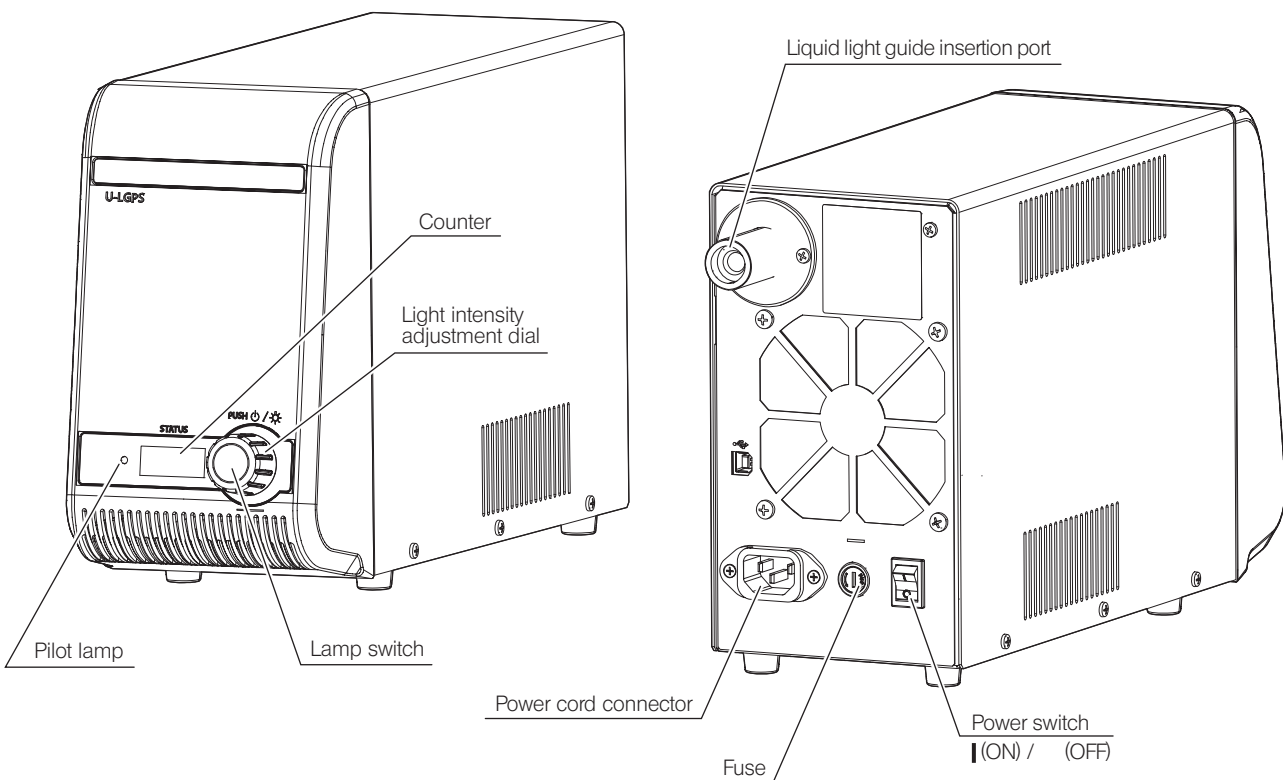

Power supply unit  
U-RFL-T

TIP


For details, see the instruction manual provided with the U-RFL-T.

Xenon lamp housing  
U-LH75XEAP0  
Power supply unit  
U-RX-T

TIP

For details, see the instruction manual provided with the U-LH75XEAP0 and U-RX-T.

Tube lens unit  
MVX-TLU

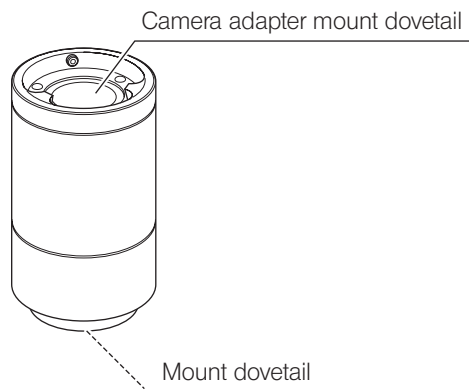



# 3 Summary of transmitted light brightfield observation procedure

## 3-1 Preparation

1. Set the shutter knob to "●" to close the shutter. (Page 15)
2. Mount the fluorescence mirror unit(s) that match the desired observation. (Page 36)
3. Turn the mirror unit turret to engage the desired fluorescence mirror units in the light path. (Page 13)
4. Turn on the mercury burner and wait until the arc stabilizes (for about 5 to 10 minutes). (Page 14)  
(If the burner is not centered, center it.)

## 3-2 Observation procedure

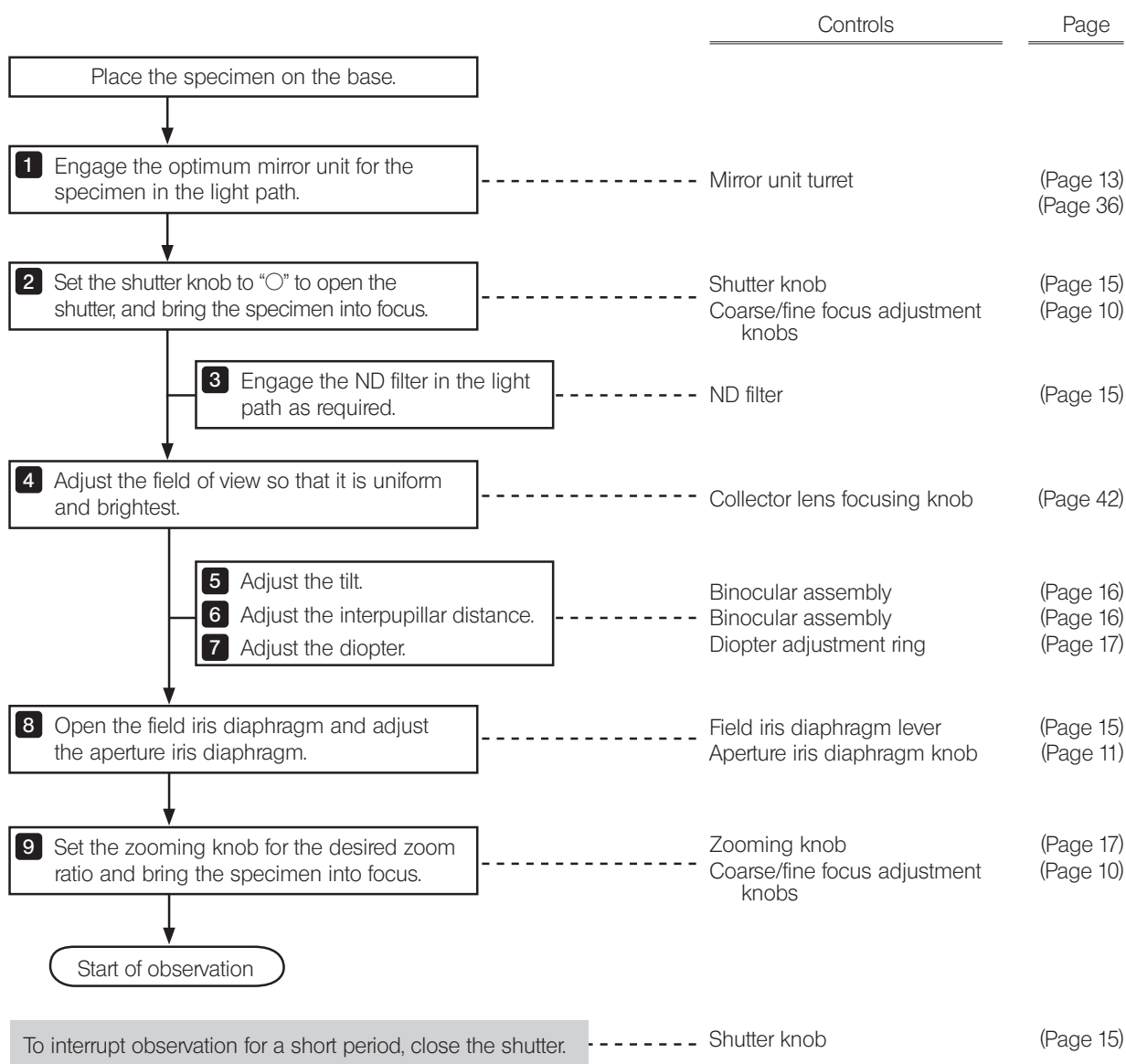

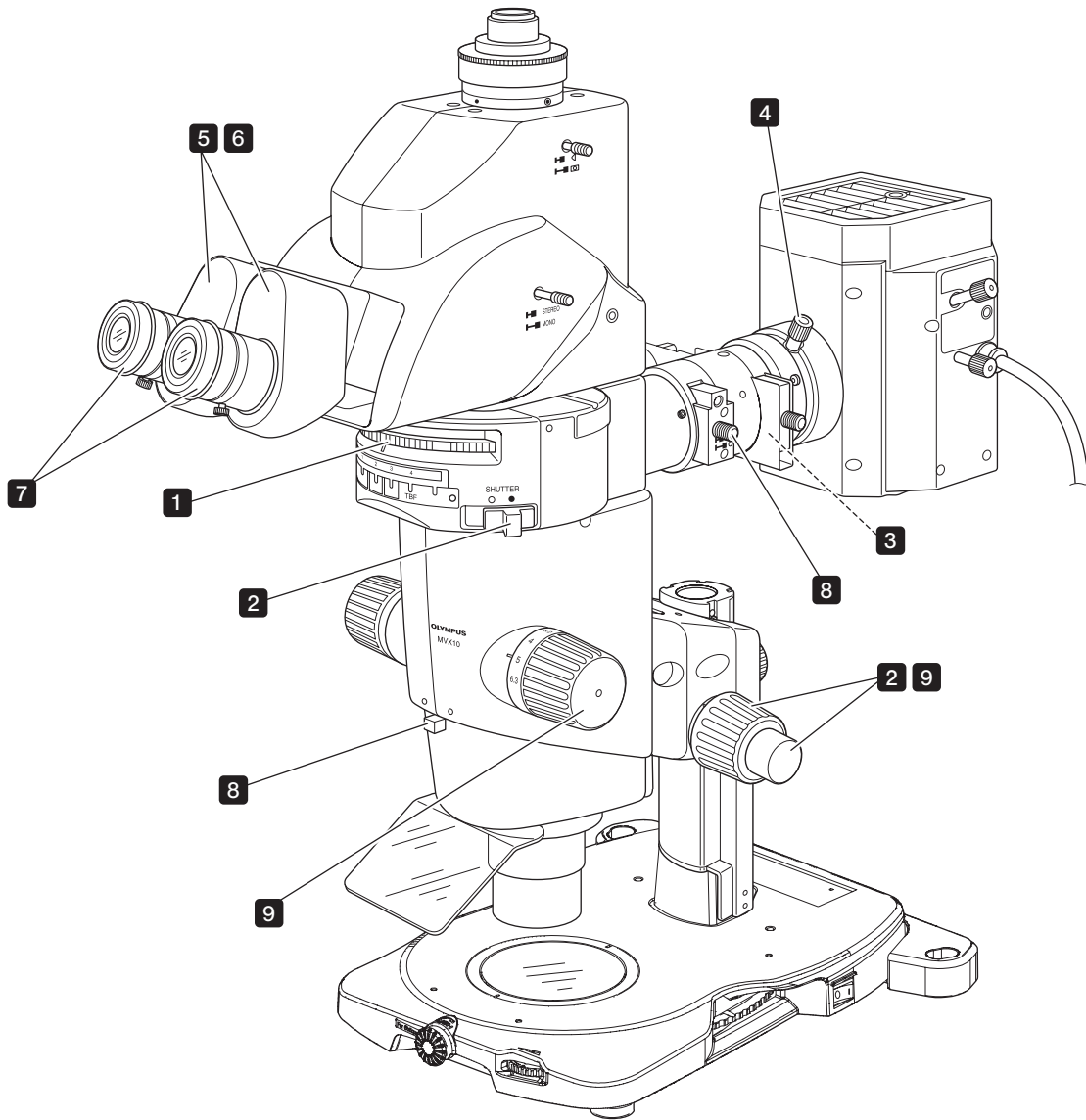

**TIP** Make a photocopy of this spread and post it near the microscope for quick reference.


For the observation by transmitted light illumination, refer to the instruction manual provided separately.

# 4 Operation

## 4-1 Base

For the operation of the base, refer to the instruction manual provided separately.

## 4-2 Microscope body and focusing assembly

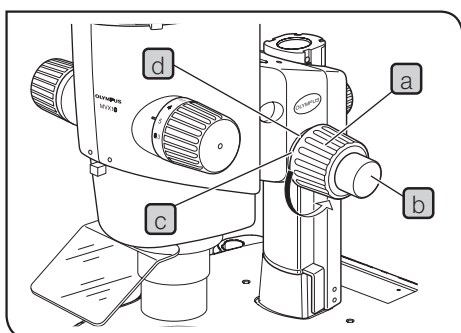

### 1 Adjusting the focus

With both the coarse focus adjustment knob **a** and fine focus adjustment knob **b**, rotating the knob in the direction of the arrow lowers the microscope body.

- Stroke per turn of the coarse focus adjustment knob: 36.8 mm
- Stroke per turn of the fine focus adjustment knob: 0.77 mm

### 2 Adjusting the rotation tension of the coarse focus adjustment knob

**NOTE** The rotation tension of the coarse focus adjustment knob can be adjusted with the rotation tension adjustment ring **c**. Do not rotate the knobs on the left and right in opposite directions, for this will damage the internal mechanism.

**TIP** This adjustment is intended to facilitate the rotation of the knob while preventing the microscope body from lowering spontaneously. For best ease of use, it is recommended to adjust the rotation tension slightly tighter than the level at which spontaneous lowering of the microscope body occurs.

**1** Rotate the rotation tension adjustment ring **c** by inserting the Allen screwdriver into the hole **d** on the ring periphery. Rotating the ring clockwise increases the rotation tension of the coarse focus adjustment knob, and rotating counterclockwise decreases it.

**NOTE**


- If the microscope body falls down by its own weight or the focus obtained by fine focusing is lost immediately, the rotation tension adjustment may be too light. In this case, rotate the ring clockwise to increase the rotation tension.
- If the rotation tension adjustment is too tight, delicate focusing will be impossible and the knob may be damaged. Particularly, never rotate the fine focus adjustment knob quickly while its rotation tension is extremely tight.

### 3

## Engaging and disengaging the zooming knob click stop position

#### TIP

- When the click stop knob is engaged to ON, the click stop function is engaged for each zoom ratio indicated with the zooming knob. When the knob is disengaged to OFF, the zoom magnification can be varied continuously and finely near the click groove.
- Click stop can be provided for the nine zoom ratios from 0.8 to 5. The click stop is engaged when the microscope leaves the factory.

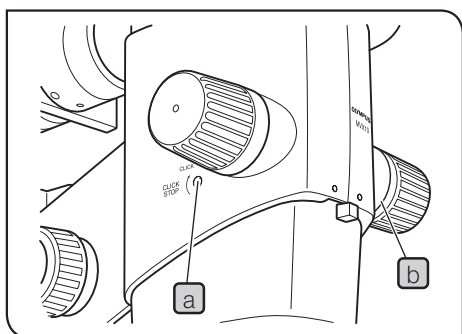

#### To disengage the click stop function

Rotate the click stop ON-OFF screw **a** counterclockwise (by about two turns from the ON position, in the direction opposite to the arrow) using the Allen screwdriver.

#### NOTE

**Do not rotate too much, or the cover may be damaged.**

#### To engage the click stop function

Rotate the click stop ON-OFF screw **a** fully clockwise (in the direction of the arrow) using the Allen screwdriver. The zooming knob then stops at every position corresponding to the zoom ratio indicated on the zoom ratio indicator **b**.

### 4

## Adjusting the aperture Iris diaphragm

#### TIP

Adjusting the aperture iris diaphragm improves the insufficiency of brightness in the peripheral section light at intermediate zoom ratios as well as the depth of focus. However, setting the aperture iris diaphragm too narrowly degrades resolution.

When performing fluorescent light observation, it is recommended to open the aperture iris diaphragm fully for bright observation.

#### 1


Adjust the aperture iris diaphragm knob **a** to the left or right.


#### TIP

Rotating the ring toward the left "O" opens the aperture; rotating it toward the right "⊗" closes it. Adjust while monitoring the observed image to confirm the contrast and focal depth improvement effects.

#### NOTE

**Do not close the aperture too much, for this may cause degradation in resolution and/or insufficiency of brightness in the peripheral section.**

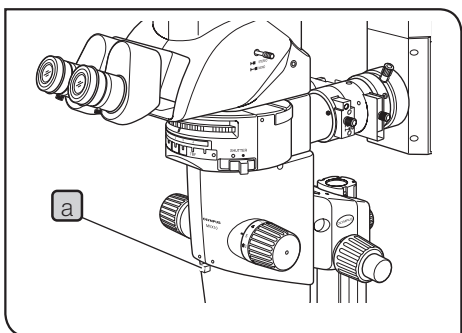

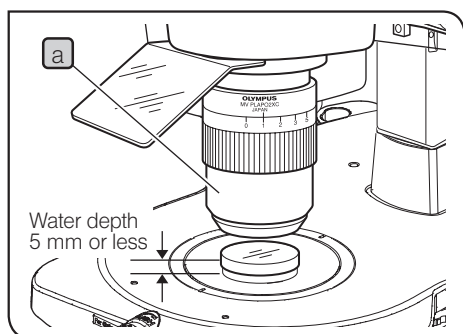

## 5 Using the objective correction collar

### TIP

- The MVPLAPO 2XC objective is equipped with a correction collar **a**, which corrects the aberration produced by a medium such as water or plastic container. When observing the specimen through a liquid or petri dish cover, rotate the correction collar to find the position that provides the best contrast.
- Correction is possible for aberrations equivalent to a water depth of about 5 mm.
- The effect of the correction collar may be less sensible when the zoom ratio is low or the aperture iris diaphragm is stopped down.
- When the MVX-2RE revolving nosepiece is used, hold the objective so that the revolving nosepiece does not displace from the position when turning the correction collar **a**.

## 4-3 Coaxial fluorescence illuminator


### 1 Selecting the fluorescence mirror unit

**TIP** Use the fluorescence mirror unit that is optimum for each purpose of observation.


Mirror unit name	Dichroic mirror	Excitation filter	Barrier filter	Application
U-MCFPHQ/XL	DM450HQ	BP425-445HQ	BA460-510HQ	ECFP
U-MGFPHQ/XL	DM485HQ	BP460-480HQ	BA495-540HQ	EGFP
U-MYFPHQ/XL	DM505HQ	BP490-500HQ	A515-560HQ	EYFP
MRFPHQ/XL	DM565HQ	BP535-555HQ	BA570-625HQ	RFP
U-MGFP/XL	DM505	BP460-490	BA510IF	EGFP (U-MGFPA/XL For dye separation)
U-MGFPA/XL			BA510-550	

#### Spectral characteristics of the filters


1. Filter for fluorescent protein CFP U-MCFPHQ/XL

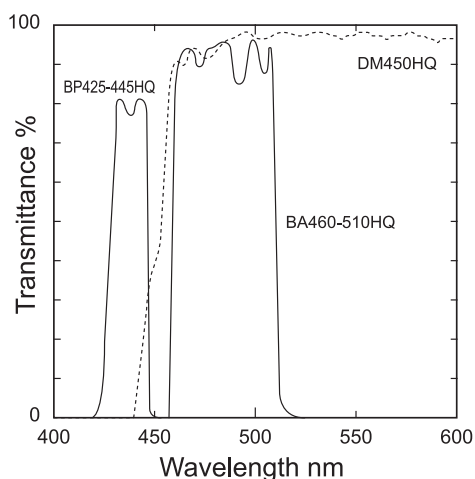


2. Filter for fluorescent protein GFP U-MGFPHQ/XL


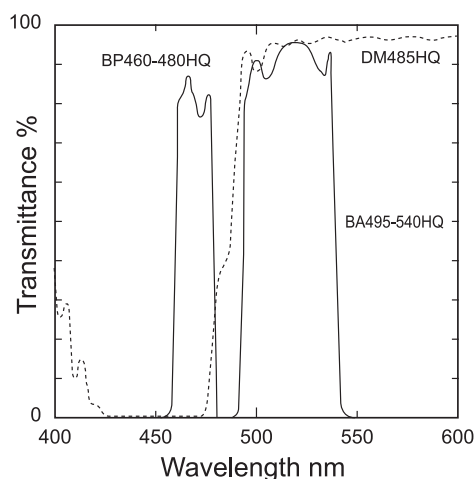


3. Filter for fluorescent protein YFP U-MYFPHQ/XL

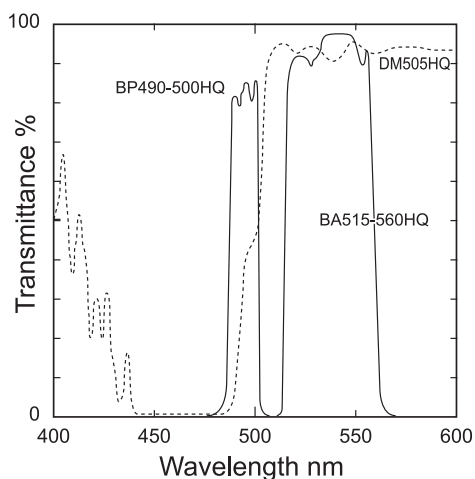

4. Filter for fluorescent protein RFP U-MRFPHQ/XL

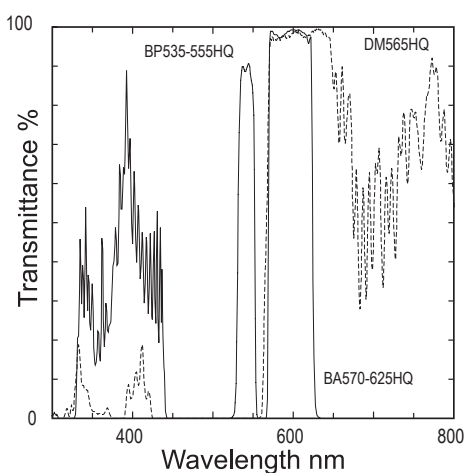

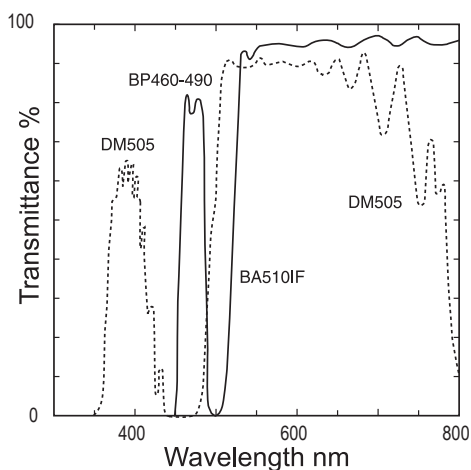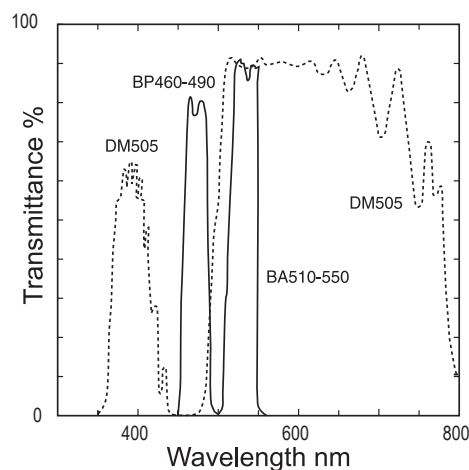5. IIB excitation filter (wide band) U-MGFP/XL6. IIB excitation filter (wide band) U-MGFPA/XL

## 2 Turning the burner ON


**NOTE** The burner should be centered before the first use after it has been mounted or replaced (see page 42). When resetting the burner hour counter, press and hold the reset button until the reading becomes "0.0".

**1** Set the main switch of the power supply unit to "I" (ON). The arc image of the burner will stabilize in 5 to 10 minutes after the burner is ignited.

**TIP** A discharge-type mercury burner may not turn on by the first try due to the characteristics of the burner. If a burner does not turn on, set the main switch to "O" (OFF), wait for 5 to 10 seconds and set the main switch to "I" (ON) again.

**NOTE**


- To avoid shortening the service life of the burner, do not turn a burner off for 2 hours after turning it on but use the shutter in this period.
- When turning on a mercury burner that has been turned off, wait for about 10 minutes after it is turned off, because it cannot be turned on unless the mercury vapor in the burner has cooled down and liquefied.
- If the lamp housing is opened while the burner is on, the power supply stops to ensure safety. In this case, set the main switch to "O" (OFF) and wait for more than 10 minutes before retrying to turn it on again. Do not open the lamp housing unless it has cooled down sufficiently.

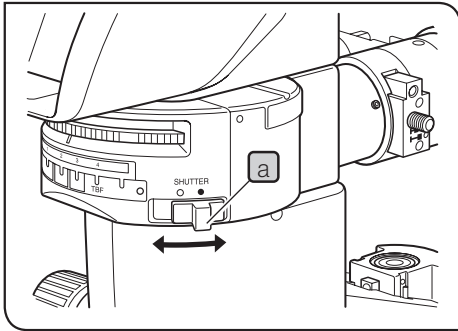

### 3 Opening/closing the shutter

**TIP** The shutter can be used to close the shutter temporarily when changing the mirror unit turret or interrupting observation for a short period.


- 1 When the shutter knob **a** is set to “●”, the shutter is engaged to shut off the excitation light.
- 2 When the shutter knob **a** is set to “○”, the shutter is disengaged and opened.

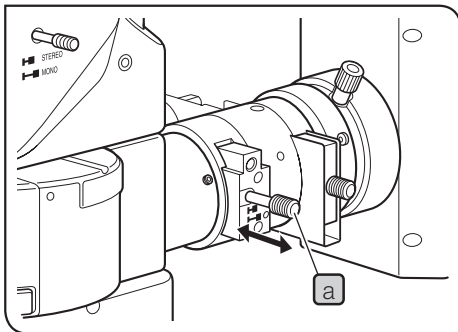

### 4 Using the field Iris diaphragm

**TIP** • The centering of the field iris diaphragm can be adjusted (see page 43).

- Push in or pull out the field iris diaphragm lever **a** as required to prevent color fading due to fluorescence in parts other than the observed region.

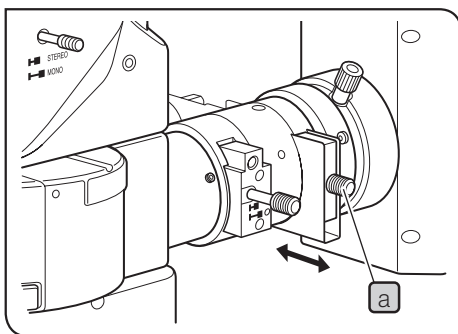

### 5 Switching the filter slider knob


**TIP** • Hold the filter slider knob **a** of coaxial fluorescence illuminator and move it toward the left or right to select one of the two filter pockets or idle position.

- The idle position is selected when the knob is pulled to the rightmost stop position. Pushing the knob at the click position one step after the idle position selects filter pocket 1, and pushing it further to the leftmost stop position selects filter pocket 2.
- To adjust the brightness using filters, mount the required ND filter (32ND6, 32ND12, 32ND25, 32ND50) or frost filter (32FR80) in the filter pockets and select the desired filter pocket with the filter slider knob.

#### **CAUTION**

When the mercury burner has been turned on for an extended period, the metallic part of the filter slider knob and the ND filter get very hot. Be careful when switching the filter slider knob or changing the filter.

## 4-4 Observation tube

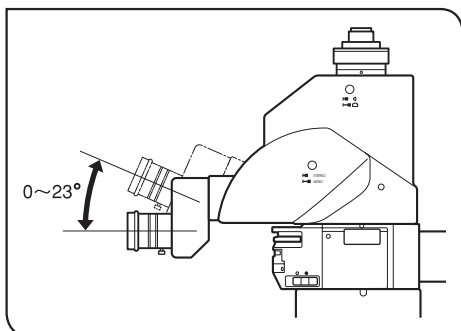


### 1 Adjusting the tilt

Adjust the height and tilt of the eyepieces of the binocular assembly to the most comfortable viewing position

- 1 Holding the binocular assembly with both hands, raise or lower it to the desired position.


**NOTE** Do not attempt to force the binocular assembly past the upper or lower stop positions. Applying excessive force could destroy the mechanism.

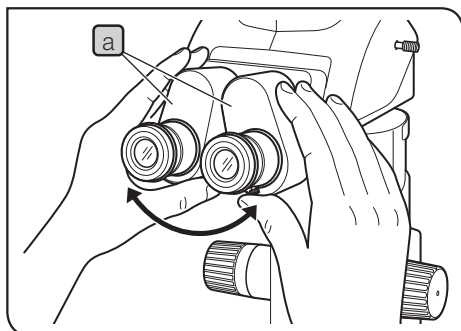

### 2 Adjusting the interpupillar distance

**NOTE** Be sure to hold the binocular assembly **a** with both hands to make this adjustment.

- 1 While looking through the eyepieces, hold the left and right of the binocular assembly **a** and adjust by opening or closing it for binocular vision until the left and right fields of view coincide completely.

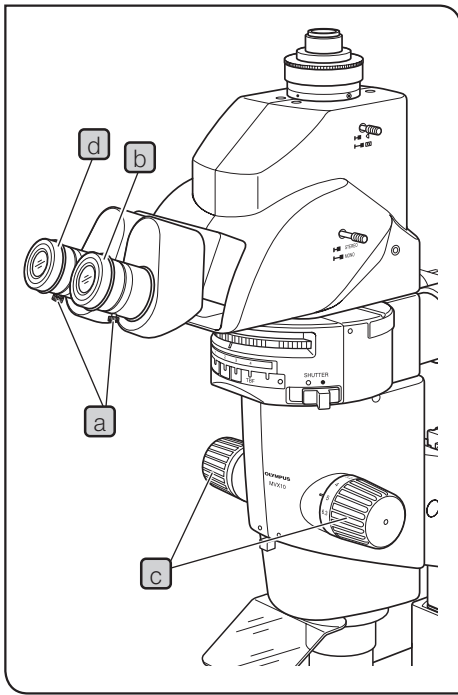

### 3 Adjusting the diopter (zoom parfocality adjustment)


**TIP** Make sure that the eyepiece clamping knobs **a** are tightened. Adjustment of the diopter according to the user's eyesight ensures parfocality even when the zoom magnification is changed.

- 1** Looking into the right eyepiece, rotate the diopter adjustment ring **b** so that the outer periphery of the field of view is as sharp as possible.
- 2** Place an easy-to-observe specimen on the stage.
- 3** Rotate the zooming knob **c** to a low zoom ratio and focus the specimen using the coarse and fine focus adjustment knobs.
- 4** Rotate the zooming knob **c** to the highest zoom ratio and focus the specimen using the coarse and fine focus adjustment knobs.
- 5** Rotate the zooming knob **c** to the lowest zoom ratio, then focus the specimen by rotating the left and right diopter adjustment rings **d** instead of the coarse and fine focus adjustment knobs.

#### When using the micrometer eyepiece

- 1** Look into the micrometer eyepiece and focus on the micrometer scale by rotating the diopter adjustment rings **b**.
- 2** Place an easy-to-observe specimen on the stage.
- 3** Rotate the zooming knob **c** to a low zoom ratio and focus on the specimen looking into the micrometer eyepiece and using the coarse and fine focus adjustment knobs.
- 4** Rotate the zooming knob **c** to the highest zoom ratio and focus on the specimen using the coarse and fine focus adjustment knobs.
- 5** Rotate the zooming knob **c** to the lowest zoom ratio, then focus the specimen by rotating only the diopter adjustment ring **d** of the nonmicrometer eyepiece instead of the coarse and fine focus adjustment knobs.

**TIP** Note (or memorize) the diopter readings of the left and right eyepiece scales so that they can be duplicated quickly in the next observation.

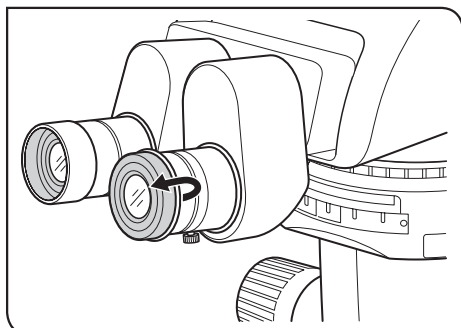


#### 4 Using the eye shades

When wearing eyeglasses

Use with eye shades in their normal folded-down position. This will prevent eyeglasses from being scratched by the eyepieces.

When not wearing eyeglasses

Extend the folded eyeshades in the direction of the arrow. This makes observation easier by preventing the inverse incidence of light from between the eyepieces and your eyes.

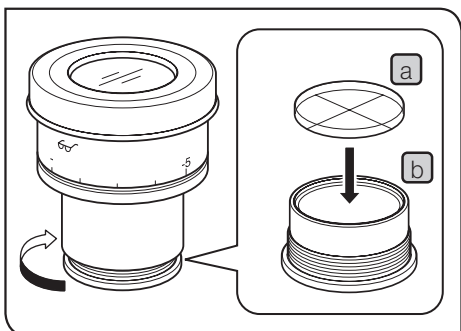


#### 5 Mounting the eyepiece micrometer disk

When the WHN10X-H eyepieces are used, an eyepiece micrometer disk **a** can be mounted. Use 24 mm dia. x 1.5 mm thick micrometer disks.

**1** Following figure, turn the built-in micrometer-mounting frame **b** to the arrow direction to remove it from the eyepiece and place a micrometer disk **a** into the mounting frame so that the surface with the model indication faces downward.



**TIP** The micrometer-mounting frame may be tight for certain micrometer disks. In this case, turn the frame by holding the circumference with a light, uniform force or by applying the frame against a rubber sheet. Do not grasp the frame with a strong force, as this may deform the frame and make it harder to remove it.

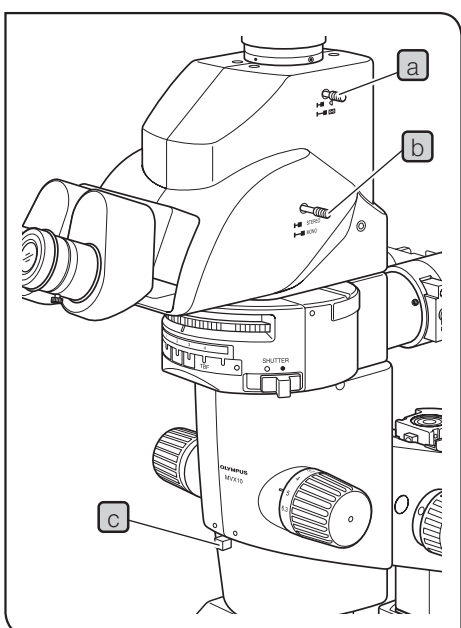**NOTE** Be careful not to touch the lens surface with your finger.



#### 6 Selecting the light path

Slide the light path selector lever **a** to select the desired light path.

Light path selector lever	Indication	Intensity Ratio
Pushed in		Binocular 100%
Pulled out		Camera 100%

#### 7 Switching the monaural/stereo view

Slide the mono/stereo selector lever **b** for stereo observation based on pupil division.

Mono/stereo selector lever	Indication	Observation
Pushed in*	 STEREO	Stereo
Pulled out	 MONO	Standard

\* The brightness of the peripheral section of the field becomes insufficient when the aperture iris diaphragm knob **c** is stopped down.

# 5 Camera observation and photomicrography

When camera observation or photomicrography is required, use the MVX-TTRS tilting trinocular head or MVX-TLU tilting lens unit.


The MVX-TV 1XB/MVX-TV 1XC/TV 0.63XC-2 exclusive camera adapter can be installed on this microscope.

When the MVX-TV 0.63XC-2 camera adapter is used, insufficiency of brightness in the peripheral section may sometimes be noticeable at low to intermediate zoom ratios. In this case, stop down the aperture iris diaphragm knob on the front of the microscope body.

## 1 Selecting the camera adapter magnification


Set the magnification of the camera adapter according to the size of the CCD in the digital camera.

The following figures show the camera observation areas when the WHN10X-H eyepieces with a field diameter of 22 are used.

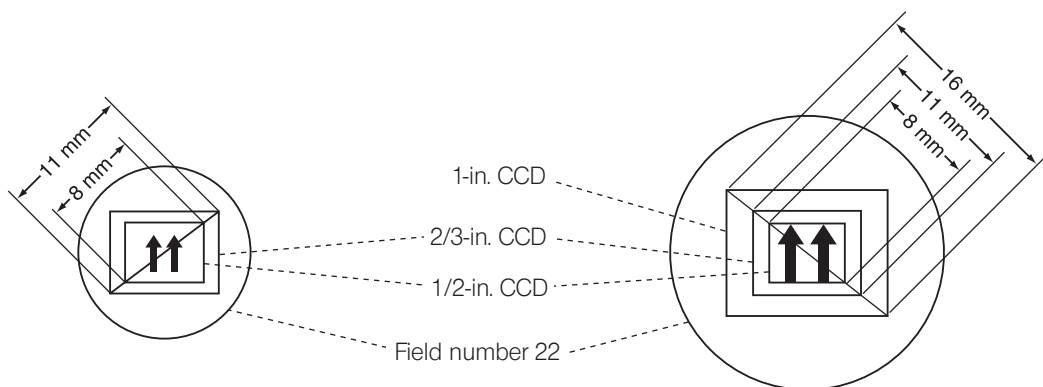

When the 0.63X camera adapter is used  
(The field may be cut off when the 1-inch CCD is used)

When the 1X camera adapter is used

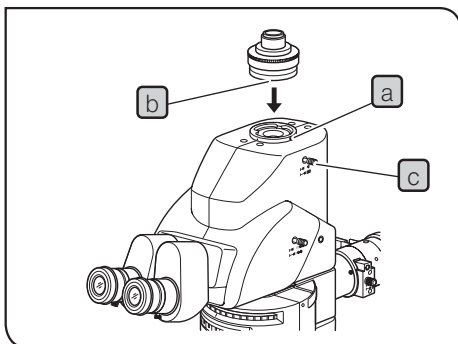

## 2 Attaching the camera adapter

- TIP**
- Attach the camera adapter to the camera first, and mount the camera adapter to the tube.
  - For procedures to attach the camera adapter to the camera, refer to Instruction Manual provided with the camera.

- 1 Using the Allen screwdriver, fully loosen the straight tube clamping screw **a** of the straight tube mount on the top of the tilting trinocular head.
- 2 Fit the round dovetail **b** of the camera adapter into the straight tube mount of the trinocular head, and tighten the clamping screw **a**.


## 3 Selecting the camera light path

Pull out the light path selector knob **c** to select the Camera 100% light path setting.

## 4 Adjusting the parfocality between observation image and monitor image


**TIP** The parfocality adjustment makes correction of focusing unnecessary when the observation image is switched to the image monitored.

- 1** Looking into the eyepieces, select a high zoom ratio and focus on the specimen.
- 2** Select the camera light path to view the monitored image and select a low zoom ratio.

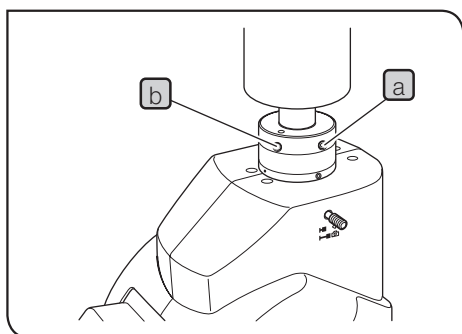


When using the MVX-TV 0.63XC-2

- 3** Using the Allen screwdriver, loosen the parfocality adjustment clamping screw (LOCK) **a** on the camera adapter.
- 4** While observing the monitored image, turn the parfocality adjustment screw (FOCUS) **b** slowly to focus on the specimen.
- 5** When the specimen is focused, tighten the clamping screw **a** using the Allen screwdriver.

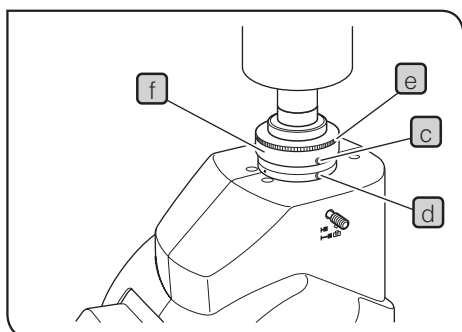

When using the MVX-TV 1XC

- 3** Using the Allen screwdriver, loosen the parfocality adjustment clamping screw **c** and camera adapter clamping screw **d**.
- 4** While observing the monitored image, hold the upper part **e** of the camera adapter and rotate the lower part **f** to focus on the specimen.
- 5** When the specimen is focused, tighten the clamping screws **c**, **d** using the Allen screwdriver.

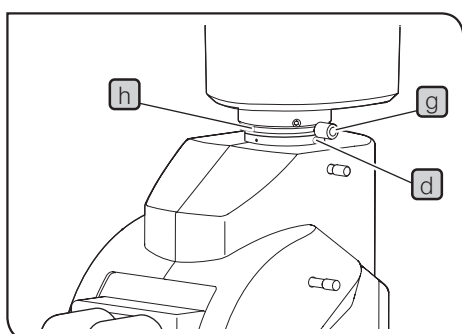

When using the MVX-TV 1XB

- 3** While observing the monitor image, adjust its focus by rotating the parfocality adjustment area **g** of the camera adapter.
- 4** When the specimen is focused, tighten the clamping screw **h**.

### Adjusting the camera angle

- 1.** Loosen only the camera adapter clamping screw **d** and rotate the camera to adjust its angle.
- 2.** After adjusting the angle, tighten the clamping screw **d**.

# 6 Troubleshooting

Under certain conditions, performance of this unit may be adversely affected by factors other than defects. If a problem occurs, please review the following list and take remedial action as needed. If you cannot solve the problem after checking the entire list, please contact your local EVIDENT representative for assistance.

Problem	Cause	Remedy	Page
1. Incomplete binocular vision.	Interpupillary distance is not correctly adjusted.	Adjust it correctly.	16
	Diopter adjustment is incomplete.	Complete diopter adjustment.	17
2. Field of view is cut off or illuminated unevenly.	Stereo observation is performed and the aperture iris diaphragm is stopped down excessively.	Open the aperture.	11
	The trinocular head and intermediate attachment are not correctly mounted.	Mount them correctly.	38, 39
	Light path selector knob is stopped midway.	Set correctly to the desired position.	18
	Mirror unit turret or filter slider knob is in an intermediate position.	Set them in the correct positions.	15
3. Dust is visible in the field of view.	Dust on the specimen.	Remove dust.	3
	Dust on eyepiece.	Remove dust.	
4. Excessive image contrast.	Aperture is stopped down excessively.	Open the aperture to proper diameter.	11
5. Resolution problems: • Image is not sharp. • Insufficient contrast.	Objective is not correctly mounted.	Mount it correctly until it is stopped.	35
	Dust on objective front lens.	Clean lens surfaces	3
	Dust on top or lower lens of zoom microscope body.		
6. Specimen image blurs when zoom ratio is changed.	Diopter of the eyepieces is not correctly adjusted.	Adjust it correctly.	17
	Not in complete focus on specimen.	Focus specimen correctly at a high zoom ratio.	17
7. Coarse focus adjustment knobs rotate with too much resistance.	Rotation tension adjustment ring is too tight or a module is attached on the microscope body.	Loosen the ring properly.	10
8. Zoom microscope body drops or specimen goes out of focus during observation.	Rotation tension adjustment ring is too loose.	Tighten the ring properly.	10
9. Image looks doubled.	The MVX-CA2X is set to an intermediate position	Push or pull the lever into a stop position.	26
10. Background noise is noticeable in the fluorescence.	The excitation light is too strong.	Insert the ND filter and set the excitation light to the minimum required intensity level.	38

# 7 Specifications

Item	Specification
1. Zoom microscope body MVX-ZB10	Zoom magnification system. Zoom drive system: Horizontal knob. Click stop ON-OFF switchable per zoom ratio. Zoom ratios: 10 (0.63X to 6.3X). Zoom ratio indications: 0.63, 0.8, 1, 1.25, 1.6, 2, 2.5, 3.2, 4, 5, 6.3. Objective mount: Threaded mount Built-in aperture iris diaphragm.
2. Objective	MVPLAPO 0.63X : NA 0.15, WD 87 mm, FN 22. MVPLAPO 1X : NA 0.25, WD 65 mm, FN 22. MVPLAPO 2XC : NA 0.5, WD 20 mm, FN 22., with correction collar.
3. Eyepiece	WHN10X-H*: FN 22, with diopter adjustment ring. * $\varnothing$ 24 mm dia., 1.5 mm thick eyepiece micrometer disk can be inserted.
4. Focusing assembly SZX2-FOFH	Focusing system: Rack & pinion roller guide (with coarse focus adjustment knob rotation tension adjustment ring). With built-in gas spring counter balance. Coaxial coarse/fine focus adjustment knobs. Focusing stroke: 80 mm. Coarse knob stroke per turn: 36.8 mm. Fine knob stroke per turn: 0.77 mm.
5. Tilting trinocular head MVX-TTRS	Tilting angle: 0° to 23° Light path selection: 2 steps (Bi 100% or Camera 100%). Interpupillary distance adjustment: 51 to 76 mm. Eyepiece clamping knob provided. Eyepiece: WHN10X-H
6. Coaxial fluorescence illuminator MVX-RFA	Illumination magnification: 1X (FN 22). Observation switching: Mirror unit turret. Number of mountable fluorescence mirror units: 3 (with idle positions). Field iris diaphragm provided (centering possible). Shutter provided. Filter slider knob provided (2 filter pockets + 1 idle position). Available observation methods: Reflected light fluorescence and transmitted light.
7. Lamp housing for mercury burner	U-LH100HG: Lamp housing for 100 W mercury burner. U-LH100HGAP0: Apo lamp housing for 100 W mercury burner. Mercury burner: USH-103OL (Ushio Inc.) or HBO103W/2 (OSRAM) Power Supply Unit: U-RFL-T
8. Operating environment	Indoor use. <ul style="list-style-type: none"> <li>• Altitude : Max. 2000 meters (6440 ft).</li> <li>• Ambient temperature: 5° to 40°C (41° to 104°F).</li> <li>• Relative humidity: 80% for temperatures up to 31°C (88°F), decreasing linearly through 70% at 34°C (93°F), 60% at 37°C (99°F) to 50% relative humidity at 40°C (104°F).</li> <li>• Supply voltage fluctuations: <math>\pm</math>10%.</li> <li>• Pollution degree: 2 (in accordance with IEC60664).</li> <li>• Installation (overvoltage) category: II (in accordance with IEC60664)</li> </ul>

# 8 Optical characteristics

Total magnification/actual field

- Total magnification = Objective power x Zoom ratio x Eyepiece magnification.

- Actual field area =  $\frac{\text{Eyepiece FN}}{\text{Objective power} \times \text{Zoom ratio}}$


Objective	Eyepieces	
	WHN10X-H (Field number: 22)	
	Total magnification	Actual field (mm)
MVPLAPO 0.63X	4X to 40X	55.4 to 5.54
MVPLAPO 1X	6.3X to 63X	34.9 to 3.49
MVPLAPO 2XC	12.5X to 125X	17.4 to 1.74

# 9 Operation of other modules


## 9-1 Revolving nosepiece MVX-2RE

The revolving nosepiece accepts two MVPLAPO series objectives to enable observation using a wider variety of magnifications. The objectives can be switched quickly by simply rotating the revolving nosepiece horizontally.

### 1 External view

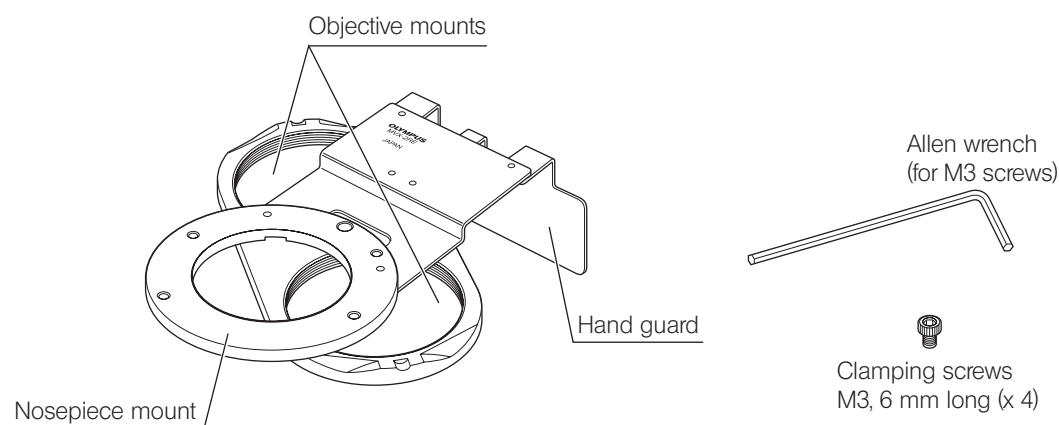


### 2 Assembly

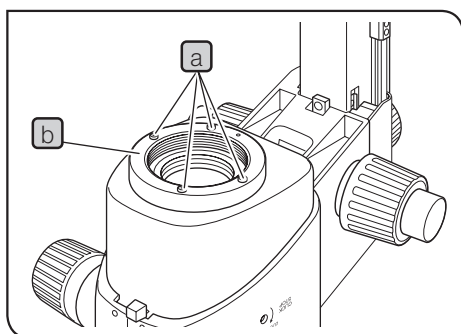

- 1 Remove the trinocular head and coaxial fluorescence illuminator from the zoom microscope body.
- 2 Remove the zoom microscope body and focusing assembly from the pillar, and place them upside down on a flat desk surface.

**NOTE** Place a soft sheet made of rubber or similar material on the desk surface.


- 3 Remove the objective from the zoom microscope body and then remove the objective mount (b) by loosening the four clamping screws (a) using the provided Allen wrench (for M3 screws).

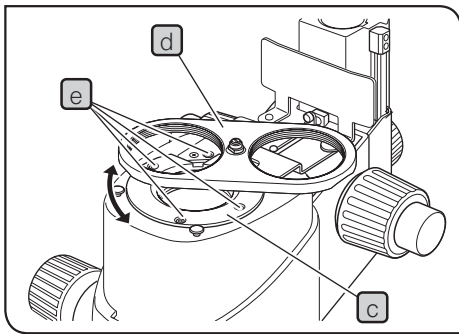

- 4** Place the nosepiece mount **c** (with its objective mounts **d** facing upward) where the objective has been by aligning the screw holes. Using the Allen wrench (for M3 screws), clamp the nosepiece mount with the provided four clamping screws (M3, 6 mm long) **e**.


**NOTE**

- As the screw holes may be hidden behind the objective mounts **d**, clamp the nosepiece mount while rotating it.
- Before installing the focusing assembly on the pillar, adjust the coarse and fine focus adjustment knobs so that the hand guard **e** does not protrude from the lower end of the focusing assembly **f**.

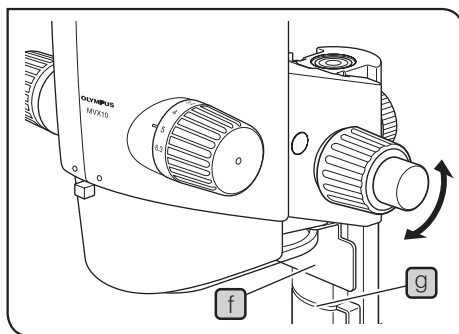

- 5** Install the focusing assembly on the pillar, and install the observation tube and coaxial fluorescence illuminator which have been removed above in the original positions.
- 6** Attach two objectives onto their respective objective mounts **d** by screwing.

### 3 Operation

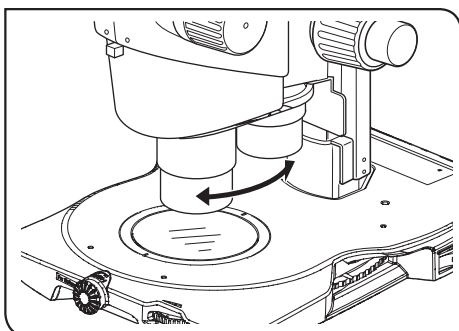

- 1** Hold the objectives and gently rotate the nosepiece until a click position where the objective to be used is engaged in the light path.

### 4 Caution

- When transporting the microscope, do not hold it by the revolving nosepiece.
- The parfocal property when the objective is switched is not completely guaranteed.

## 9-2 Magnification changer MVX-CA2X

When the magnification changer has been installed on the MVX-TTRS tilting trinocular head by the dealer, the observation magnification can be switched between 1X and 2X with a selector lever.

Pushed-in position: 1X


Pulled-out position: 2X

Be sure to push in or pull out the selector lever in a position where it is stopped securely.

**NOTE** Not to spoil the optical performance, always have your dealer perform the assembly work.

### Assembly (operation performed by the dealer)

**NOTE** The MVX-CA2X is shipped with grease applied on it. Be careful not to stain the lenses with the grease during assembly.

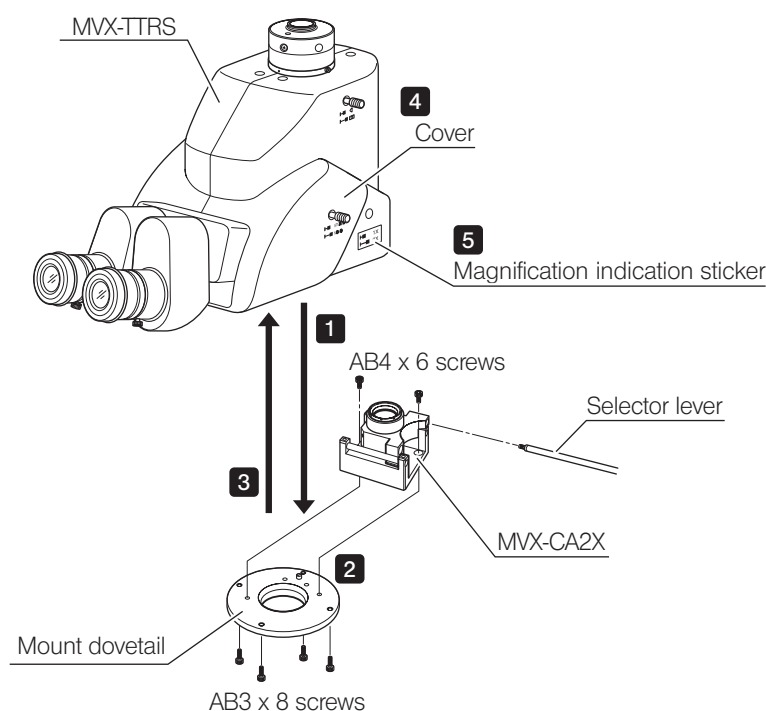

- 1** Using the Allen wrench, remove the four AB3 x 8 screws from the mount dovetail on the MVX-TTRS.
- 2** Using the two AB4 x 6 screws and the Allen screwdriver, attach the MVX-CA2X onto the mount dovetail by aligning the positioning pin.
- 3** Place the mount dovetail with the MVX-CA2X on the original position by aligning the positioning pin, and then clamp using the four AB3 x 8 screws.
- 4** Remove the cover of the lever insertion hole of the MVX-TTRS and screw the selector lever into it.
- 5** Attach the magnification indication sticker below the lever.


## 9-3 BX stage adapter type 1 SZX-STAD1

This adapter is for installation of a U-SRG or U-SRP rotary stage on the SZX2-STL large base or the SZX2 illumination base. When the U-SRP rotary stage is used together with the U-FMP mechanical stage, X-Y directional movement becomes possible, which is convenient for framing during photomicrography and camera observation.

### 1 Applicable bases & restrictions

Base	Applicable objectives	Restrictions
<ul style="list-style-type: none"> <li>• Large base SZX2-STL</li> </ul>		None.
<ul style="list-style-type: none"> <li>• Quad position LED transmitted light illumination base SZX2-ILLTQ</li> <li>• Single position LED transmitted light illumination base SZX2-ILLTS</li> </ul>	MVPLAPO 0.63X MVPLAPO 1X MVPLAPO 2XC	STAD mount for ILLT SZX2-STADM is necessary for using the stage adapter. (Refer to the instruction manual for the SZX2-ILLTQ/ILLTS LED transmitted light illumination base.)


### 2 Assembly

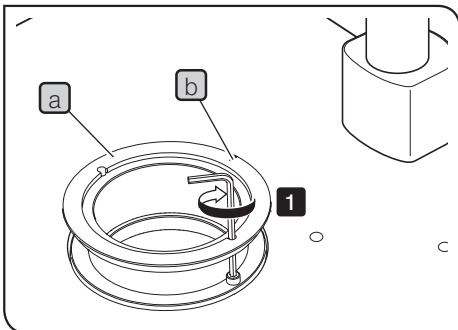

**TIP** Use the tools shown in the table below.

Tool	Description
Allen wrench (for M4 screw)	Provided with SZX-STAD1
Hexagon socket screw (M4)	Provided with SZX-STAD1

- Place the notch **b** of SZX-STAD1 **a** on the back side of the base, and secure the screws (2 pcs.) using the Allen wrench to attach SZX-STAD1 **E** to a base.

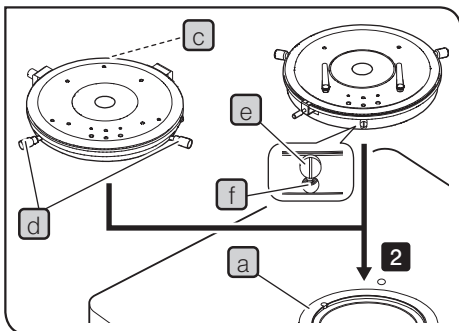

- Attach U-SRP or U-SRG2 to SZX-STAD1 **a**.

#### Attaching U-SRP

Place the positioning pin (cylinder type) **c** on the back side of the base, and rotate the centering knob **d** in clockwise direction to secure it.

#### Attaching U-SRG2


Place the protrusion **e** on the front side of the base, and rotate the clamping screw **f** in clockwise direction to secure U-SRG2 using the allen screwdriver provided with the microscope frame.

## 9-4 Stage adapter type 1 SZH-STAD1

This adapter has the same function as the SZX-STAD1 BX stage adapter type 1, but the usable stages with this adapter is only the BH2-SH horizontal knob stage.

### 1 Assembly


**TIP** STAD mount for ILLT SZX2-STADM is necessary for using the SZX2-ILLTQ/ILLTS LED transmitted light illumination base. For details, refer to the instruction manual for the SZX2-ILLTQ/ILLTS LED.

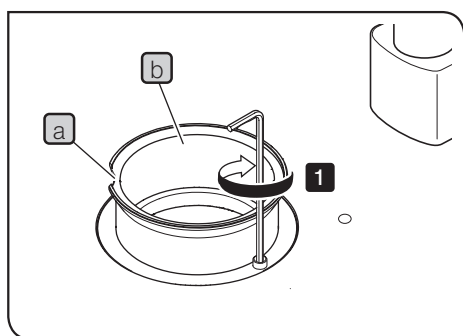

**TIP** Use the tools shown in the table below.

Tool	Description
Allen wrench (for M4 screw)	Provided with SZH-STAD1
Hexagon socket screw (M4)	Provided with SZH-STAD1

**1** Placing the notch **a** on the front side of the base, secure the screws (2 pcs.) using the Allen wrench to attach SZX-STAD1 **b** to a base.

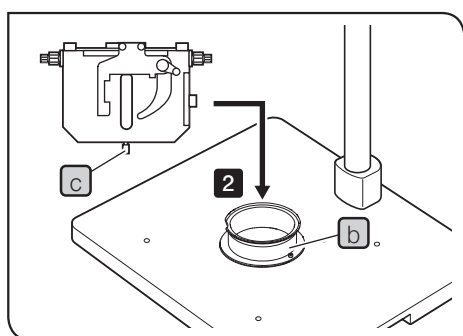

**2** Place the stage clamping knob **c** of BH2-SH (horizontal knob stage) on the front of the base, rotate the clamping knob **c** in clockwise direction to secure it, and attach BH2-SH.

## 9-5 BX stage adapter type 2 SZX-STAD2


This adapter is for installing the U-SIC4 large stage\* on the SZX2-STL large base\*\*. When this stage adapter is used, the provided SZH-P400 auxiliary pillar should be used to cover the height of the stage adapter.

\* The U-SVL or U-SVR BX stage can also be used but the operability is degraded, and the U-SVLB and U-SVRB cannot be used due to the long stage knobs.


\*\* STAD mount for ILLT SZX2-STADM is necessary for using the SZX2-ILLTQ/ILLTS LED transmitted light illumination base. In addition, the LED transmitted illumination built in SZX2-ILLTQ/ILLTS cannot be used for observation. For details, refer to the instruction manual for the SZX2-ILLTQ/ILLTS LED.

### 1 Assembly

**NOTE** For simplified transmitted light observation, place the mirror knob on the front and use a frosted filter.

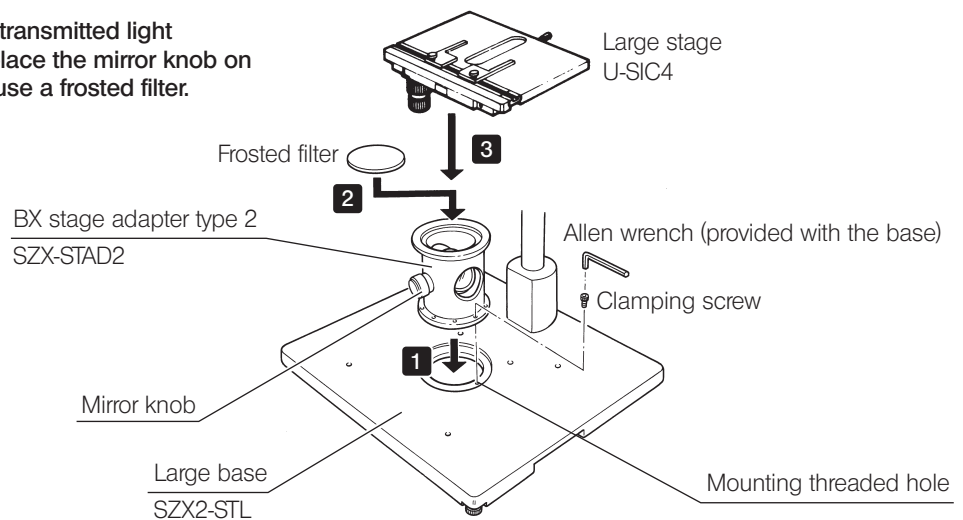

### 2 Simplified transmitted light observation

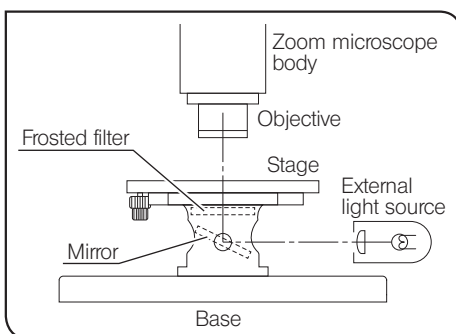

- 1** Illuminate the specimen with an external light source (LSD illuminator, light guide illuminator, etc.). Light the external light source as shown in the figure on the left and irradiate the mirror assembly.
- 2** Eliminate irregularities in illumination.
  - 1) Align the microscope body center with the center of the SZX-STAD2 adapter.
  - 2) Set the zooming knob on the microscope body to the minimum magnification and focus the stage top surface.
  - 3) While observing through the eyepiece, rotate the mirror knob to adjust the mirror angle so that the entire field of view is illuminated uniformly.


**TIP** When observing using an oblique illuminator, remove the frosted filter and, while observing through the eyepiece, tilt the mirror gradually until optimum contrast can be obtained.

### 3 Caution

- Do not project the image of the external light source filament on the frosted surface of the frosted filter. Otherwise, the frosted filter may deteriorate.
- Use neutral detergent to clean the frosted filter.
- In transmitted light observation at a total magnification of no more than 10X, the field of view may be obscured in the peripheral sections depending on the stage in use.

## 9-6 Gliding stage SZH-SG


### 1 External view & nomenclature

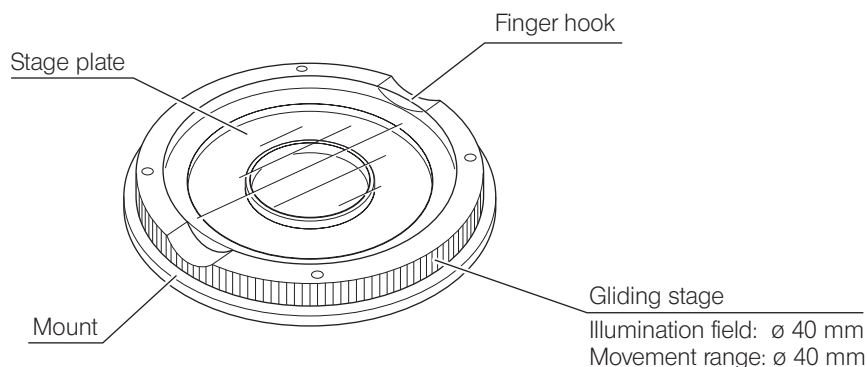

### 2 Assembly

- NOTE**
- If dirt or metallic powder is attached on the gliding surfaces, be sure to clean them.
  - When placing the gliding stage parts, be sure to avoid bringing the gliding surfaces in contact with the desk surface.

**TIP** STAD mount for ILLT SZX2-STADM is necessary for using the SZX2-ILLTQ/ILLTS LED transmitted light illumination base. For details, refer to the instruction manual for the SZX2-ILLTQ/ILLTS LED.

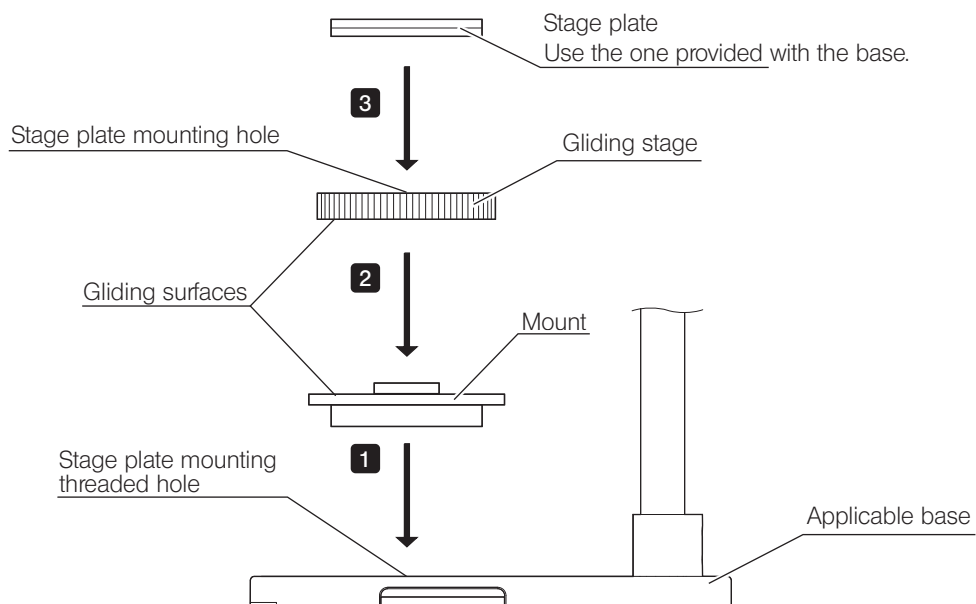

**TIP** Clean the gliding surfaces periodically.


### 3 Operation

Hold the gliding stage by the surrounding edge and move it horizontally.

## 9-7 Cup stage SZH-SC

### 1 External view & nomenclature


**NOTE** This stage is applicable only to reflected light illumination.

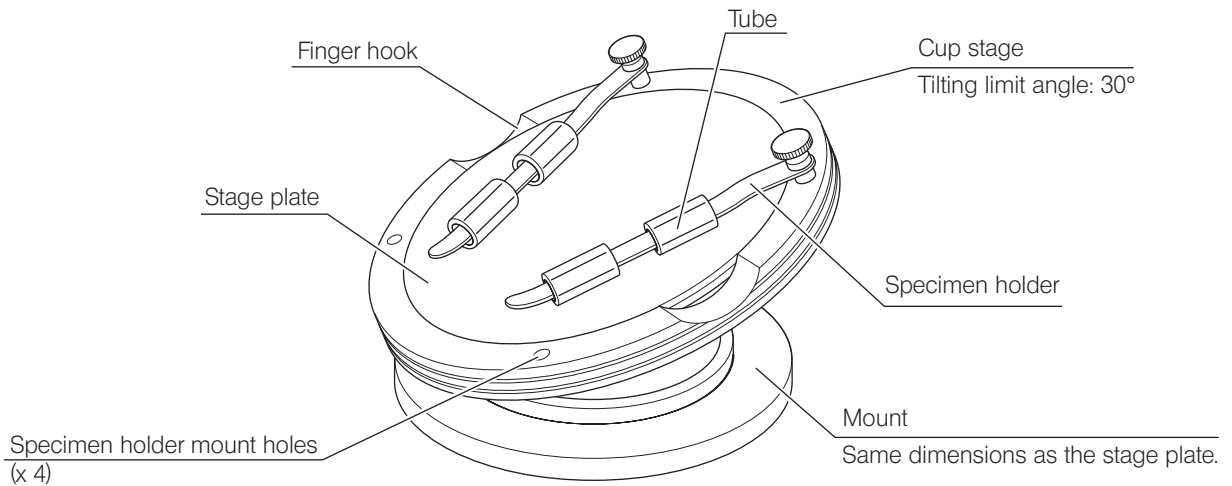

### 2 Assembly

**NOTE** Before assembly, remove dirt and dust from the parts and perform the operation cautiously.

**TIP** STAD mount for ILLT SZX2-STADM is necessary for using the SZX2-ILLTQ/ILLTS LED transmitted light illumination base. For details, refer to the instruction manual for the SZX2-ILLTQ/ILLTS LED.

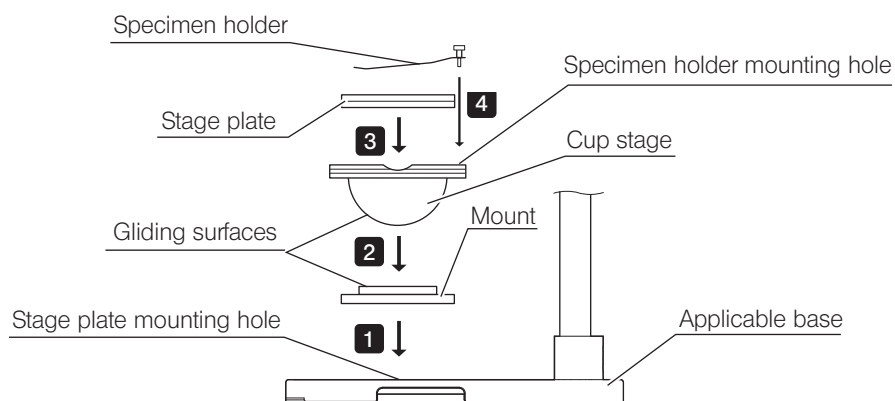


- 1 Fit the cut stage mount into the stage plate mounting hole of an applicable base.
- 2 Place the cup stage on the mount.  
Before placing, clean the gliding surfaces on the cup stage and mount by wiping with a clean cloth.

3 Attach the stage plate.

4 Attach the specimen holder.


**TIP** Clean the gliding surfaces periodically.

### 3 Operation

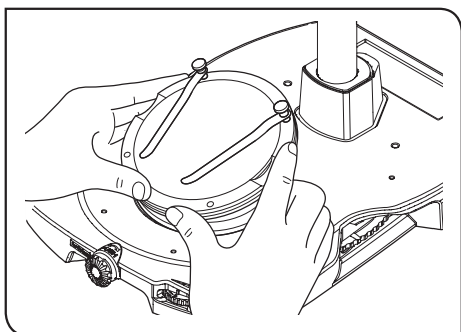

- 1** Place a specimen on the stage plate, hold the edge of the cut stage and tilt it slowly.

**TIP** If the specimen slips off, fix it using the provided specimen holder.

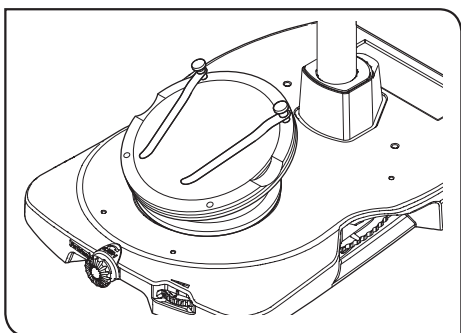

**TIP** To fix a container such as a petri dish, fit the provided tube around the specimen holder and hold the container with it.

- NOTE**
- Take care not to touch the gliding surfaces of the cup stage and mount with a hand. If they are contaminated, clean with neutral detergent before use.
  - The cup stage may move spontaneously if a load of 20 g or more is applied to its outer periphery.
  - When the cup stage is tilted while a specimen with large height is placed on it, the image may be defocused. If this happens, adjust focusing again.


# 10 Assembly

## 10-1 Assembly diagram

The diagram below shows how to assemble the various modules. The numbers indicate the order of assembly.


**NOTE**

**When assembling the microscope, make sure that all parts are free of dust and dirt, and avoid scratching any parts or touching glass surfaces.**

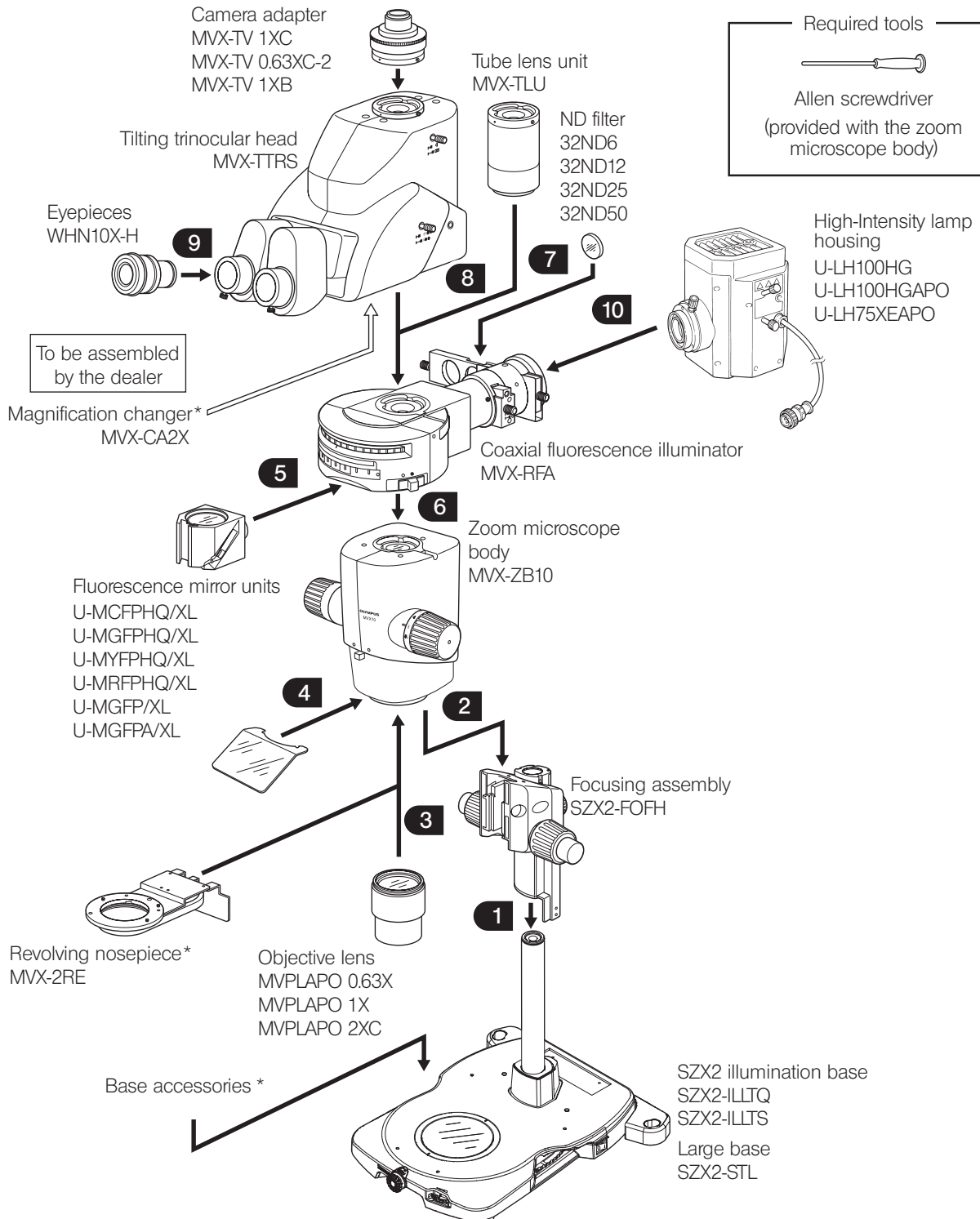

\* The assembly methods are described in "9 Operation of other modules" on page 24.

## 10-2 Detailed assembly procedures

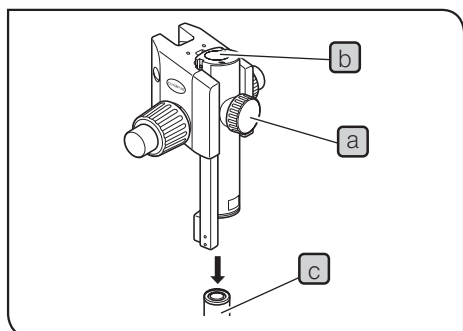


### 1 Mounting the focusing assembly

- 1 First loosen the focusing assembly clamping knob **a** completely and, while holding the focusing assembly with both hands, insert the pillar **c** into the mounting hole **b** from below.


**NOTE** Insert slowly. Do not apply excessive force.

- 2 Lower the focusing assembly until it stops, then tighten the focusing assembly clamping knob **a**.

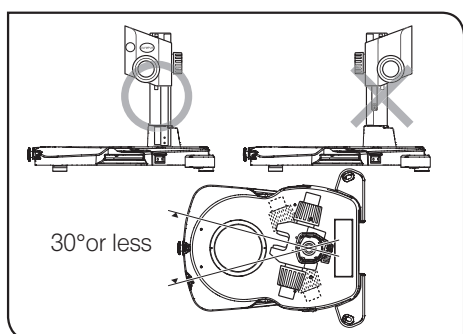

**CAUTION** To prevent the microscope from turning over, the focusing assembly must be installed as shown in the illustration marked "O" in figure, and its pivot angle must be limited to 30°. If the focusing assembly is placed on the wrong side, the microscope will turn over.

**NOTE** If the clamping knob **a** is tightened while the pillar **c** is not completely inserted into the mounting hole **b**, the plate spring supporting the pillar from the inside will deform and the pillar will not be able to penetrate into the hole.

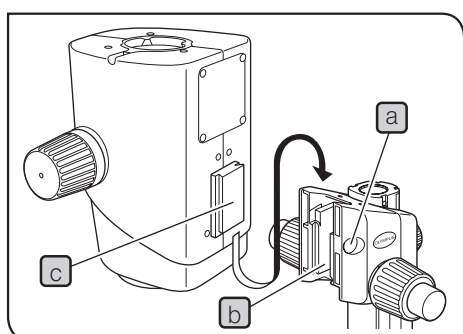


### 2 Mounting the zoom microscope body

- 1 Remove the cap **a** on the focusing assembly by inserting a thin object into the notch.
- 2 Using the provided Allen screwdriver, loosen the dovetail mount clamping screw inside the cap on the focusing assembly by rotating it by 2 or 3 turns (counterclockwise).
- 3 Gently insert the dovetail mount **c** on the rear of the microscope body into the dovetail mount **b** on the focusing assembly.

**NOTE** Do not insert the mount at an angle or with excessive force, for this may cause malfunctions.

- 4 When the microscope body has been inserted until it stops, tighten the clamping screw using the Allen screwdriver.
- 5 Place the cap **a** in the original position.

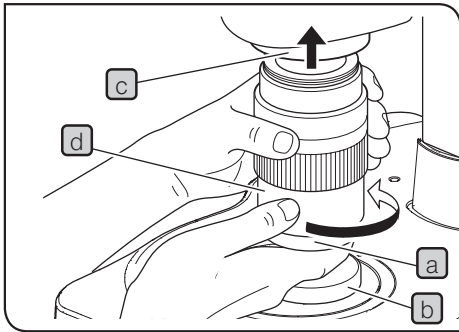


### 3 Mounting the objective

**NOTE** Since the objective is heavy, there is a risk of dropping it during mounting or dismounting. To prevent this, take the following countermeasures.



- Attach the provided cap **a** on the objective.
- Protect the base against the dropped objective, place the other cap **b**, a notebook or a protecting material on the base.

- 1 Mount the objective **d** on the objective mount thread **c** by rotating the objective in the direction of the arrow.

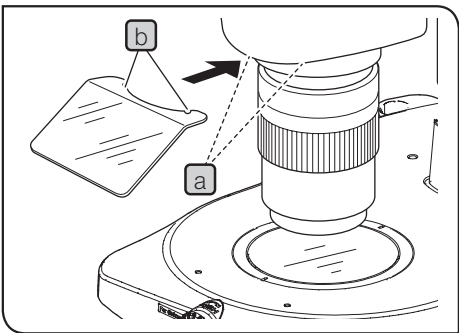

### 4 Attaching the UV shield plate

- 1 Align the mounting groove **b** of the UV shield plate with the two mounting pins **a** on the lower part of the zoom microscope body and fit the groove with a strong force.

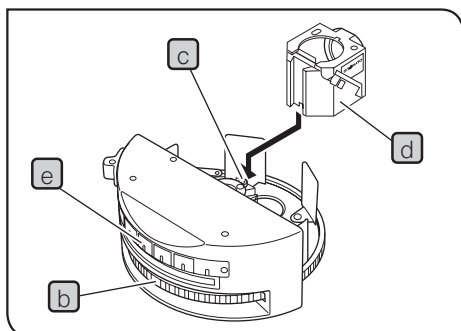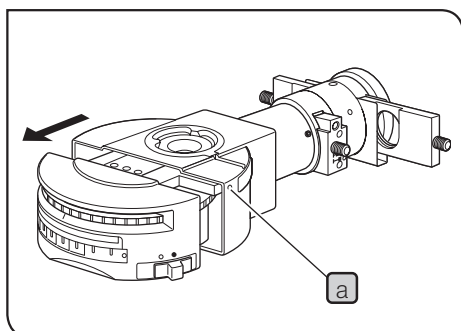

## 5 Mounting the fluorescence mirror units

Fluorescence mirror units can be mounted by in every other dovetail, or in positions with turret numbers 1, 2 and 3, of the six dovetails of the mirror unit.

**1** Using the Allen screwdriver, loosen the clamping screw hole **a** on the right end of the illuminator.

**2** Slide out the turret and place it upside down.

**TIP** After completion of assembly of the microscope, it is difficult to slide out the turret because the trinocular head comes in the way. In this case, hold the two sides of the turret with two hands and pull.

**3** Rotate the mirror unit so that the mirror unit position number (one of 1, 2 or 3) comes at the turret number indicator **b**.

**4** Loosen the clamping screw **c** using the Allen wrench.

**5** Insert the desired mirror unit **d** all the way into the dovetail position so that the abbreviation of the mirror unit name is upside down, and then tighten the clamping screw **c** firmly,

**NOTE** If the clamping screw **c** is loose, it may interfere with the inner surface of the turret cover, making impossible to rotate the turret.

**6** Select the optimum indicator plate from the provided indicator plate sheet, and insert the plate upside down in the mirror unit indicator pocket **e**.

**TIP** If the optimum indicator plate is not present, write the original name in a blank indicator plate using an oil-ink pen.

**7** Mount the other required mirror units by repeating the above steps.

**8** Place the turret in the original position and tighten the clamping screw **a** while pushing the turret.

### How to fabricate an optional mirror unit


A custom ordered or commercially available mirror unit can be fabricated by assembling a barrier filter, excitation filter and dichroic mirror, all of which are available commercially, in the U-MF/XL mirror unit frame.

However, the barrier filter and dichroic mirror should be the EVIDENT products. If a commercially available product is used, EVIDENT cannot guarantee the safety of operation.

#### Dimensional conditions of the optical parts

- Barrier filter:  $\varnothing 32_{-0.2}^{-0.1}$  mm max. thickness 4 mm
- Excitation filter:  $\varnothing 25_{-0.2}^{-0.1}$  mm max. thickness 6 mm
- Dichroic mirror: See figure on the right.

\* The arrow inscribed on the side of an EVIDENT barrier filter or excitation filter indicates the filter mounting orientation.

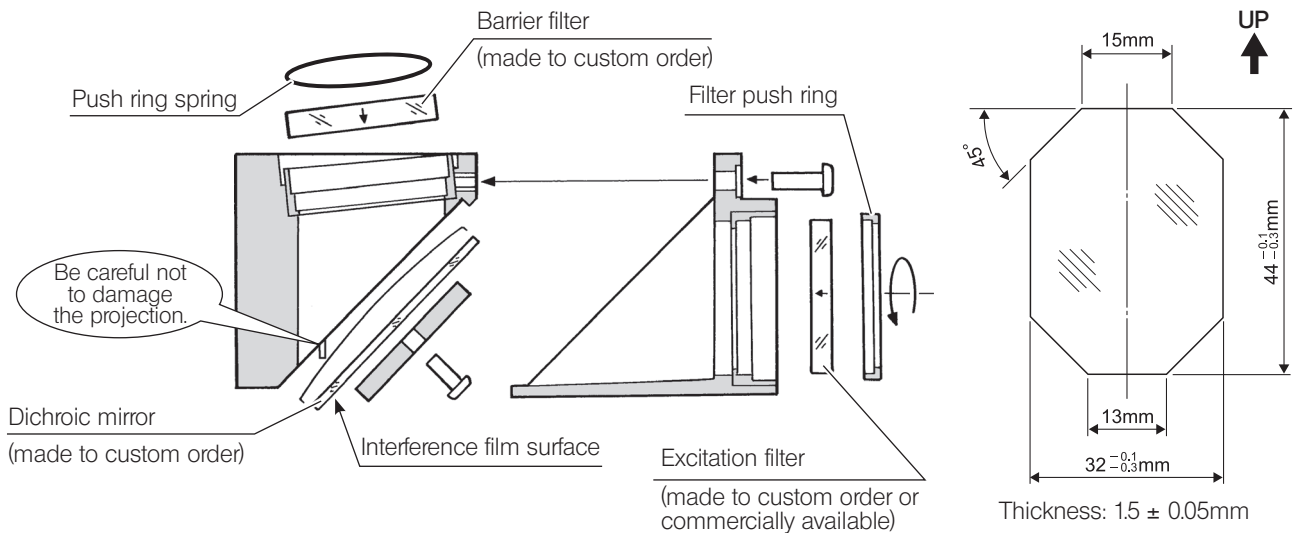


**NOTE**

When replacing the dichroic mirror, be particularly careful not to contaminate it with fingerprints, etc.

**TIP**


A mirror unit of the UIS series can also be used. However, in this case, the NA is limited in the low to medium zoom ratios due to the diameter of the filter.

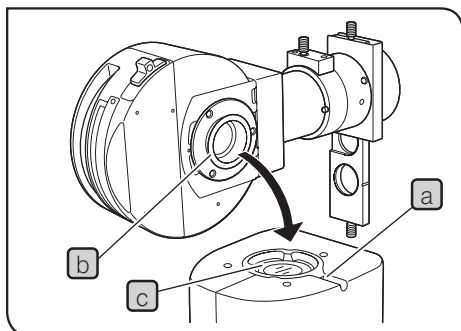

## 6 Mounting the coaxial fluorescence illuminator

- 1 Using the Allen screwdriver, loosen the illuminator clamping screw **a**.
- 2 Fit the round dovetail **b** at the bottom of the illuminator into the mount dovetail **c** on the microscope body, and tighten the clamping screw **a**.

**NOTE** To prevent the microscope from turning over after the illuminator is installed, the rotation angle should be limited to within  $\pm 60^\circ$  from the rear direction.

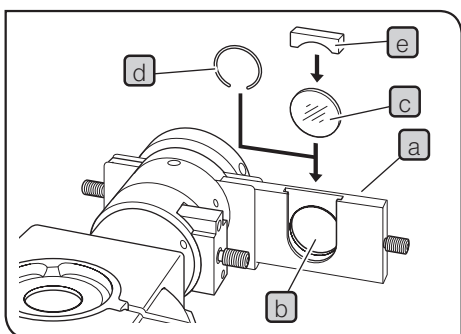

## 7 Mounting the ND filter

- TIP**
- Color fading of specimen can be delayed by attenuating the excitation light intensity using an ND filter. Use the ND filters as far as they do not hinder observation.
  - When it is difficult to mount the ND filter, detach the knob on the slider and remove the slider. Then try to mount the ND filter again.

- 1 Slide out the filter slider **a** of the coaxial fluorescence illuminator so that the filter pocket **b** is visible.

- 2 Insert the ND filter **c** in the filter pocket **b**, and insert it so that the notch of the filter push ring **d** comes on the bottom side.

**TIP** When the filter is thick (i.e. 5 mm or more), the filter will not lay down so it is not necessary to use the push ring **d**.

- 3 Apply the filter push ring **d** against the filter so that it will not lay down.

- 4 Drop the light leak prevention cap **e** into the filter slider.

**NOTE**

- Even when no filter is used, the filter push ring **d** should be mounted and the light leak prevention cap **e** should be dropped in to prevent light leak.


- Insert the filter so that the side with the indication faces the observing side. Otherwise, the filter may crack due to heat.
- Before replacing the filter, ensure that it is cooled down.

### **CAUTION**

When the mercury burner has been turned on for an extended period, the filter and its metallic parts get very hot. Be careful not to burn yourself. Also do not leave the filter slider knob in a position other than the click stop positions.


**TIP** A filter other than the ND filter can also be mounted provided that its diameter is no more than  $\varnothing 32$  mm and the thickness is no more than 6 mm.

**NOTE** Always use reflection type ND filters (32ND6, 12, 25, 50). If an absorption-type 32LND filter series or a commercially available ND filter is used, the filter crack or burn may result.

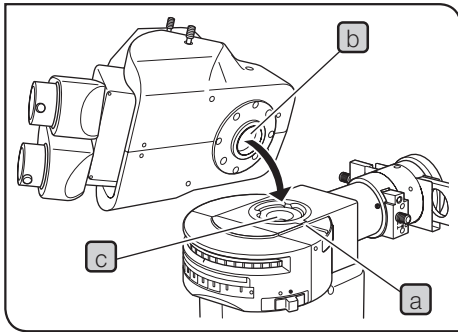

## 8 Mounting the tilting trinocular head

- 1 Using the Allen screwdriver, loosen the clamping screw **a** of the coaxial fluorescence illuminator.
- 2 Fit the round dovetail **b** on the bottom of the trinocular head into the mount dovetail **c** of the illuminator and tighten the clamping screw.

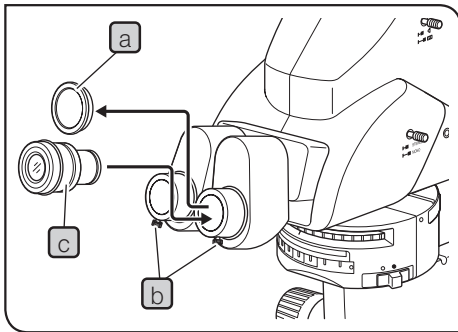


## 9 Mounting the eyepieces

- 1 Remove the eyepiece dust caps **a** and loosen the eyepiece clamping knobs **b** completely.
- 2 Gently insert the eyepiece (WHN10X-H) **c** into the left and right eyepiece sleeves all the way until they stop.
- 3 Tighten both eyepiece clamping knobs **b**.

## 10 Mounting the lamp housing for mercury burner

**TIP** For the lamp housing for xenon burner, refer to the instruction manual provided with the lamp housing. The centering method is identical to that for the mercury burner.

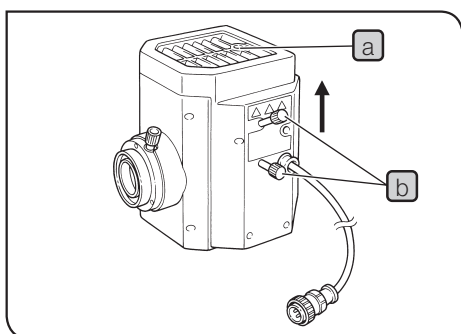

### Attaching the mercury burner

- 1 Using the Allen screwdriver, loosen the socket clamping screw **a**.
- 2 Hold the upper section of the lamp housing and pull it upward to remove the socket section.


**NOTE** To prevent malfunction, do not hold the lamp housing by the centering knobs **b**.

- 3 Place the socket section upside down as shown in figure.

**TIP** The lamp housing is equipped with the holder for transportation in the factory shipment condition or with an old burner after the burner is replaced. Remove the holder or old burner by loosening the two burner holding screws **c**.

- 4 Attach the + (positive) pole of a specified mercury burner **d** to the fixed mount on the upper side, then the - (negative) pole to the mount on the lower side.

**NOTE** Be sure to use the USH-103OL (Ushio Inc.) or the HBO103W/2 (OSRAM) mercury burner.

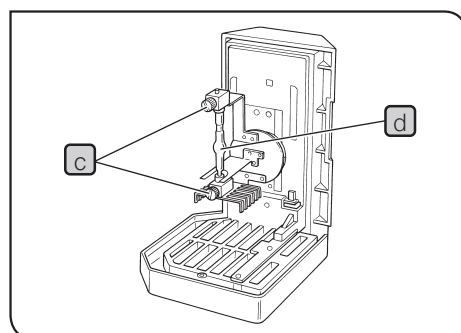

**CAUTION** Be careful and avoid leaving fingerprints or contaminants on the mercury burner. Otherwise, there is a danger of explosion due to distortion of glass caused by the stains. If the burner is contaminated, clean it by wiping gently with gauze slightly moistened with absolute alcohol.


- 5 Attach the socket section with burner to the original position and tighten the socket clamping screw **a**.  
Align the external edges of the lamp housing with those on the socket section, and push the lamp housing straight downward.

### Burner service life

USH-103OL: 300 hours

This value assumes light cycles composed of 2 hours of lighting and 30 minutes of extinction. Do not turn it on and off at a shorter cycle than the above, for this will shorten the service life of the burner.

**NOTE** After replacing the burner, reset the hour counter to "0.0" as outlined above.

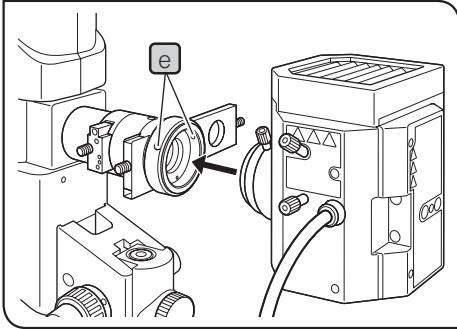




#### Attaching the lamp housing

- 1** Using the Allen screwdriver, loosen the two clamping screws **e** on the lamp housing mount hole.
- 2** Fit the lamp housing all the way as shown in figure.
- 3** Tighten the clamping screws **e** using the Allen screwdriver.

**NOTE** The surroundings of the lamp housing become very hot during and after operation. When installing the microscope system, ensure sufficient space around the lamp housing, particularly above it. Also do not mount the lamp housing obliquely.

## 10-3 Centration of the mercury (xenon) burner and field Iris diaphragm

For centering of the xenon burner, use the same procedure as the mercury burner.

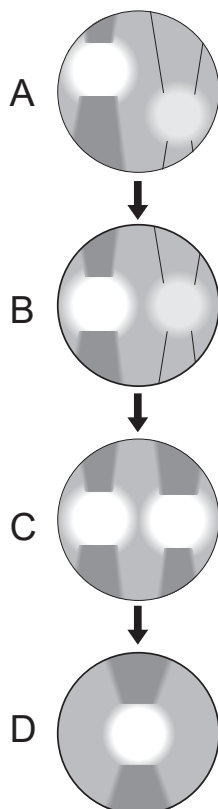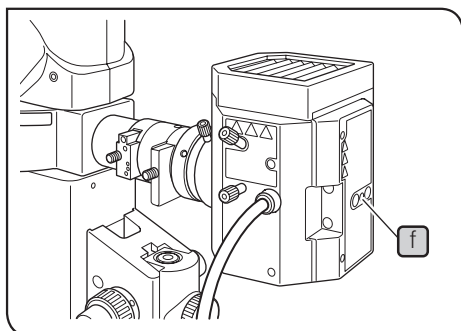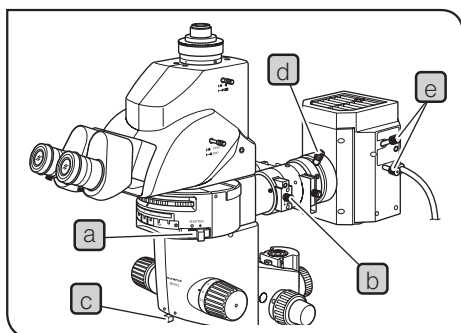

### 1 Centering the mercury burner

**TIP** Set the main switch to "I" (ON) and wait until the arc image stabilizes (for 5 to 10 minutes after ignition) before proceeding to the centering.


- 1 Set the shutter knob **a** to "●" to close the shutter.
- 2 Set the mirror unit turret to "TBF", place a sheet of white paper on the specimen surface, focus on the white sheet and set a low zoom ratio.
- 3 Engage the fluorescence mirror unit in the light path.

**CAUTION** The fluorescence illumination may contain wavelengths harmful to your eyes. The specimen must always be observed through the UV shield plate.

- 4 Remove the objective.
- 5 Pull out the field iris diaphragm lever **b** (to stop down the diaphragm).
- 6 Rotate the aperture iris diaphragm knob **c** toward the left (to open the diaphragm).
- 7 Set the shutter knob **a** to "○" to open the shutter.
- 8 Rotate the collector lens focusing knob **d** to project the arc image on the white sheet. **[A]**
- 9 Rotate the burner centering knobs **e** to bring the arc image on the center of the left (or right) half of the field. **[B]**
- 10 Insert the Allen screwdriver into the mirror focusing screw hole **f** on the rear of the lamp housing, and turn the screw to focus on the mirror-reflected image. **[C]**
- 11 Rotate the burner centering knob **e** until the arc image and mirrorreflected image overlap. **[D]**

**TIP**

- After starting observation, rotate the collector lens focusing knob **d** as required to render the observation field uniform.
- The mercury burner needs not be centered again until the time it is replaced.

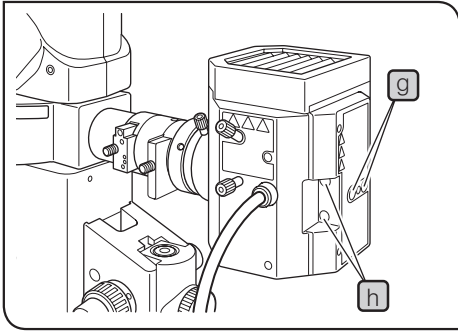


#### Precise centering of the mirror

TIP


The position of the mirror has been adjusted and locked before shipment. Only if you want more precise adjustment of the mirror position, proceed to the following steps immediately after the procedure in the previous paragraphs. Note that, once the following steps are completed, it is no longer possible to restore the mirror position in the factory shipped condition.

- 1 Using a pair of tweezers, etc., peel off the two blind stickers **g** on the rear of the lamp housing.
- 2 Fit the Allen screwdriver into each of the screws hidden below the stickers and loosen them. Loosening the two screws releases the locking of the mirror.
- 3 Peel off other two blind stickers **h** to expose the mirror centering holes.
- 4 Insert the Allen screwdriver into the screw in each mirror centering hole and adjust the centering of the mirror-reflected image.

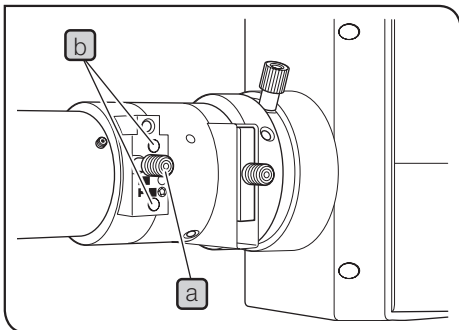

## 2 Centering the field iris diaphragm

- 1 Set the shutter knob to "●" to close the light path with the shutter.
- 2 Rotate the mirror unit turret to engage a mirror unit in the light path.
- 3 Set the shutter knob to "○" to open the shutter.
- 4 Place a specimen on the center of the base and bring the specimen into approximate focus.
- 5 Pull out the field iris diaphragm lever **a** on the illuminator to stop down the field iris diameter near the minimum size.
- 6 Insert the Allen screwdriver, provided with the microscope body, into the two field iris diaphragm centering screw **b** and adjust so that the field iris image comes on the center of the field.
- 7 Push in the field iris diaphragm lever **a** until the field iris image inscribes the field of view. If the field iris image is found to be eccentric, adjust centering again.
- 8 Open the iris until the field iris image size becomes equivalent to the field size (i.e. until the field iris image circumscribes the field of view).

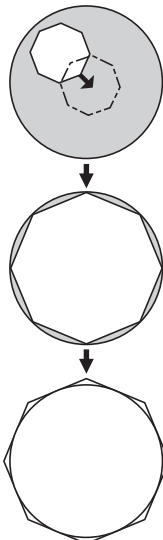

## ■ Proper selection of the power supply cord

If no power supply cord is provided, please select the proper power supply cord for the equipment by referring to "Specifications" and "Certified Cord" below:




















**Caution : In case you use a non-approved power supply cord for EVIDENT products, EVIDENT can no longer warrant the electrical safety of the equipment.**

### Specifications

Voltage rating	125 V AC (for 100-120 V AC area) or, 250 V AC (for 220-240 V AC area)
Current rating	6 A minimum
Temperature rating	60 °C minimum
Length	3.05 m maximum
Fittings configuration	Grounding type attachment plug cap. Opposite terminates in molded-on IEC configuration appliance coupling.

**Table 1 Certified cord**

A power supply cord should be certified by one of the agencies listed in Table 1 , or comprised of cordage marked with an agency marking per Table 1 or marked per Table 2. The fittings are to be marked with at least one of the agencies listed in Table 1. In case you are unable to buy locally the power supply cord which is approved by one of the agencies mentioned in Table 1, please use replacements approved by any other equivalent and authorized agencies in your country.

Country	Agency	Certification mark	Country	Agency	Certification mark
Argentina	IRAM		Italy	IMQ	
Australia	SAA		Japan	JET	
Austria	ÖVE		Netherlands	KEMA	
Belgium	CEBEC		Norway	NEMKO	
Canada	CSA		Spain	AEE	
Denmark	DEMKO		Sweden	SEMKO	
Finland	FEI		Switzerland	SEV	
France	UTE		United Kingdom	ASTA BSI	
Germany	VDE		USA	UL	
Ireland	NSAI				

**Table 2 HAR flexible cord**

Approval organizations and cordage harmonization marking methods

Approval organization	Printed or embossed harmonization marking (May be located on jacket or insulation of internal wiring)		Alternative marking utilizing black-red-yellow thread (Length of color section in mm)		
			Black	Red	Yellow
Comite Electrotechnique Belge (CEBEC)	CEBEC	<HAR>	10	30	10
Verband Deutscher Elektrotechniker (VDE) e.V. Prüfstelle	<VDE>	<HAR>	30	10	10
Union Technique de l'Electricite' (UTE)	USE	<HAR>	30	10	30
Instituto Italiano del Marchio di Qualita' (IMQ)	IEMMEQU	<HAR>	10	30	50
British Approvals Service for Electric Cables (BASEC)	BASEC	<HAR>	10	10	30
N.V. KEMA	KEMA-KEUR	<HAR>	10	30	30
SEMKO AB Svenska Elektriska Materielkontrollanstalter	SEMKO	<HAR>	10	10	50
Österreichischer Verband für Elektrotechnik (ÖVE)	<ÖVE>	<HAR>	30	10	50
Danmarks Elektriske Materialkontroll (DEMKO)	<DEMKO>	<HAR>	30	10	30
National Standards Authority of Ireland (NSAI)	<NSAI>	<HAR>	30	30	50
Norges Elektriske Materielkontroll (NEMKO)	NEMKO	<HAR>	10	10	70
Asociacion Electrotecnica Y Electronica Espanola (AEE)	<UNED>	<HAR>	30	10	70
Hellenic Organization for Standardization (ELOT)	ELOT	<HAR>	30	30	70
Instituto Portages da Qualidade (IPQ)	np	<HAR>	10	10	90
Schweizerischer Elektro Technischer Verein (SEV)	SEV	<HAR>	10	30	90
Elektriska Inspektoratet	SETI	<HAR>	10	30	90

Underwriters Laboratories Inc. (UL)  
Canadian Standards Association (CSA)

SV, SVT, SJ or SJT, 3 X 18AWG  
SV, SVT, SJ or SJT, 3 X 18AWG

# 11 Request to issue the Safety Confirmation

**NOTE** To ensure operational safety of this product, the Safety Confirmation shown on the next page must be issued by the user when the user has this product maintained, repaired or serviced by Evident Corporation.

Since Evident is responsible for ensuring the safety of our distributors, maintenance personnel and personnel, we appreciate your cooperation in issuing the Safety Confirmation.

- The user must issue the Safety Confirmation each time this product is to be maintained, repaired or serviced by Evident.
- Copy the Safety Confirmation shown on the next page, fill it out, and submit it to Evident maintenance personnel.
- If the Safety Confirmation is not issued, we may not be able to provide maintenance, repairs or other services.
- If this product is contaminated, the user shall carry out decontamination work to the fullest extent possible prior to the arrival of our maintenance personnel.

## Safety Confirmation when requesting repair/inspection

This safety confirmation is for asking you to fill out to ensure the health and safety of service personnel who pick up, repair, and/or inspect the products. Your cooperation would be highly appreciated.

Please note if the safety of service personnel cannot be ensured, your requests will not be accepted.

Please check the applicable checkboxes and circle the applicable items shown in parentheses.

Date of request	
Request details	<input type="checkbox"/> Repair ( Off-site / On-site ) <input type="checkbox"/> Inspection ( Off-site / On-site )
Product name	
Serial Number	
Biosafety level and Confirmation of disinfection/sterilization of product	<input type="checkbox"/> BSL1 <input type="checkbox"/> BSL2 <input type="checkbox"/> BSL3 <input type="checkbox"/> BSL4* <input type="checkbox"/> Not applicable <input type="checkbox"/> With my signature, I hereby confirm that this product has been disinfected/sterilized according to the disinfection/sterilization rules of our facility. <input type="checkbox"/> Disinfection/sterilization of this product has not been carried out. *We do not accept repairs/inspections of products used in the BSL4 environment.
Facility name	
Department	
Signature	
Contact details Telephone, E-mail, etc.	

Manufactured by

**Evident Corporation**

6666 Inatomi, Tatsuno-machi, Kamiina-gun, Nagano 399-0495, Japan

Distributed by

**Evident Europe GmbH**

Caffamacherreihe 8-10, 20355 Hamburg, Germany

**Evident Europe GmbH – UK Branch**

Part 2nd Floor Part A, Endeavour House, Coopers End Road, Stansted CM24 1AL, UK

**Evident Scientific, Inc.**

48 Woerd Ave, Waltham, MA 02453, USA

**Evident Scientific Singapore PTE. LTD.**

#04-04/05, 25 Ubi Rd 4, UBIX Singapore 408621

**Evident Australia PTY LTD**

Level 4, 97 Waterloo Road Macquarie Park NSW 2113, Australia

**Life science solutions**

Service Center



<https://www.olympus-lifescience.com/support/service/>

Official website



<https://www.olympus-lifescience.com>

**Industrial solutions**

Service Center



<https://www.olympus-ims.com/service-and-support/service-centers/>

Official website



<https://www.olympus-ims.com>